	<b>EN</b> Introduction <b>DE</b> Einleitung <b>ES</b> Introducción	<b>FR</b> Introduction <b>ZH</b> 简介 <b>JA</b> はじめに..... 4
	<b>EN</b> Safety Information <b>DE</b> Sicherheitshinweise <b>ES</b> Información de seguridad	<b>FR</b> Informations de sécurité <b>ZH</b> 安全信息 <b>JA</b> 安全情報 ..... 7
	<b>EN</b> Scope of delivery <b>DE</b> Lieferumfang <b>ES</b> Contenido de la entrega	<b>FR</b> Contenu de la livraison <b>ZH</b> 交货清单 <b>JA</b> 標準付属品 ..... 9
	<b>EN</b> Commissioning <b>DE</b> Inbetriebnahme <b>ES</b> Puesta en marcha	<b>FR</b> Mise en service <b>ZH</b> 调试 <b>JA</b> 試運転 ..... 10
	<b>EN</b> Pipette Check <b>DE</b> Pipettentest <b>ES</b> Comprobación de pipetas	<b>FR</b> Vérification des pipettes <b>ZH</b> 移液器检查 <b>JA</b> ピペット点検 ..... 13
	<b>EN</b> Maintenance <b>DE</b> Wartung <b>ES</b> Mantenimiento	<b>FR</b> Maintenance <b>ZH</b> 维护 <b>JA</b> メンテナンス ..... 16
	<b>EN</b> Function test <b>DE</b> Funktionstest <b>ES</b> Prueba de funcionamiento	<b>FR</b> Test de fonctionnalité <b>ZH</b> 功能测试 <b>JA</b> 機能テスト ..... 17
	<b>EN</b> Accessories and Spare Parts <b>DE</b> Zubehör und Ersatzteile <b>ES</b> Accesorios y piezas de	<b>FR</b> Accessoires et pièces de <b>ZH</b> 配件及备件 <b>JA</b> アクセサリおよびスペアパーツ ..... 20
	<b>EN</b> Disposal <b>DE</b> Entsorgung <b>ES</b> Eliminación de residuos	<b>FR</b> Mise au rebut <b>ZH</b> 处置 <b>JA</b> 廃棄 ..... 22
	<b>EN</b> Technical data <b>DE</b> Technische Daten <b>ES</b> Características técnicas	<b>FR</b> Caractéristiques techniques <b>ZH</b> 技术资料 <b>JA</b> 技術データ ..... 23
	<b>EN</b> Dimensions <b>DE</b> Abmessungen <b>ES</b> Dimensiones	<b>FR</b> Dimensions <b>ZH</b> 外形尺寸 <b>JA</b> 寸法 ..... 24
	<b>EN</b> Compliance information <b>DE</b> Informationen zur Konformität <b>ES</b> Información sobre conformidad	<b>FR</b> Information concernant la conformité <b>ZH</b> 合规性信息 <b>JA</b> コンプライアンス情報 ..... 25
	<b>EN</b> Warranty <b>DE</b> Gewährleistung <b>ES</b> Garantía	<b>FR</b> Garantie <b>ZH</b> 保修 <b>JA</b> 保証 ..... 26

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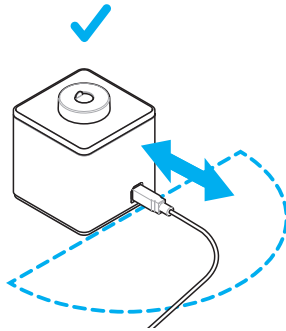
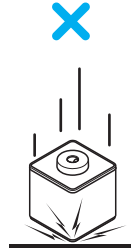
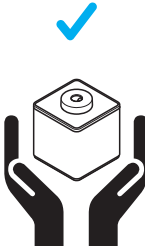
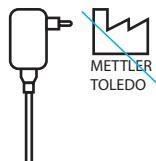
?

## EN Intended use

SmartCheck™ is a pipette check instrument. It may be used for single channel or multichannel pipettes. The intended use is to check if a pipette's delivered volume meets the set instrument tolerances. SmartCheck™ shall be used for pipette checks only and cannot be used for pipette calibration. It is only to be used with distilled water. Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

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National approval documents, e.g., the FCC Supplier Declaration of Conformity, are available online and/or included in the packaging.

► <http://www.mt.com/ComplianceSearch>

Contact METTLER TOLEDO for questions about the country-specific compliance of your instrument.

### **United States of America**

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:

1. This device may not cause harmful interference.
2. This device must accept any interference received, including interference that may cause undesired operation.

Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

The FCC Supplier Declaration of Conformity is available online.

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FR Garantie  
ZH 保修  
JA 保証

**EN Limited warranty and limitation of liability**

Seller warrants the merchandise to conform to factory specifications. As buyer's exclusive remedy in the event of a breach of warranty, seller shall repair or replace, at its option, any nonconforming merchandise or parts there of for a period on one (1) year after delivery. For more info, visit [www.mt.com/rainin](http://www.mt.com/rainin)

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**GPP®**

Good Pipetting Practice™

GPP® is a global pipette standard, ensuring consistent accuracy of pipetting processes, applicable to any pipette.

It helps to:

- Choose the appropriate pipette and tips
- Gain the range of pipetting skills necessary to produce reliable data
- Recognize the risk associated with out-of-calibration pipettes and the role of routine pipette checks vs. professional pipette service

 [www.mt.com/GPP](http://www.mt.com/GPP)

[www.mt.com/rainin](http://www.mt.com/rainin)

For more information

**Mettler-Toledo GmbH**

Im Langacher 44  
8606 Greifensee, Switzerland  
[www.mt.com/contact](http://www.mt.com/contact)

Subject to technical changes.  
© Mettler-Toledo GmbH 03/2020  
30491817A, en, de, es, fr, zh, ja

30491817