


Density Kit

for Advanced and Standard Balances

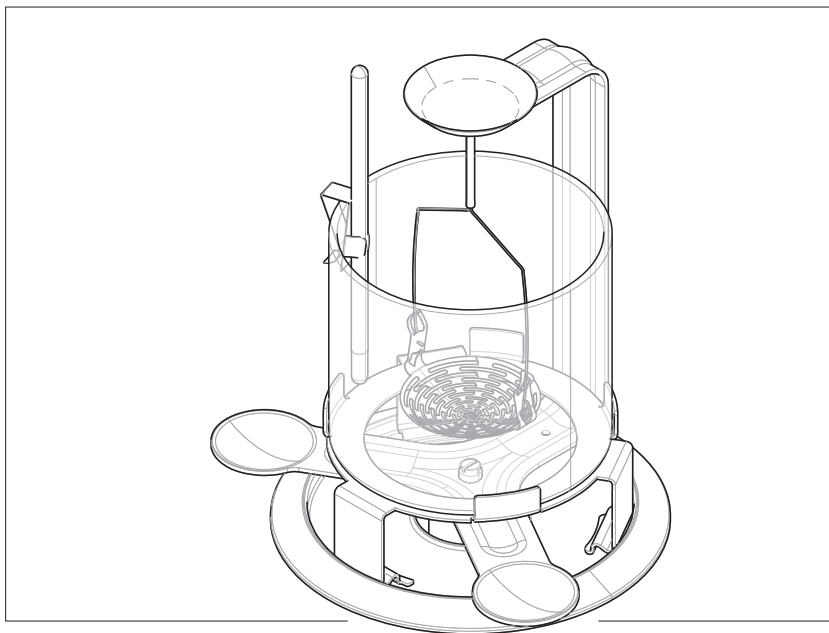

METTLER TOLEDO

Table of Contents

1	Introduction	3
1.1	Further documents and information	3
1.2	Explanation of conventions and symbols used	3
1.3	Acronyms and abbreviations	3
2	Safety information	4
2.1	Definitions of signal words and warning symbols	4
2.2	Product-specific safety notes	4
3	Installation and Putting into Operation	5
3.1	Scope of delivery	5
3.2	Compatibility table	6
3.3	Installing the density kit	7
4	Operation	8
4.1	Principle of density determination	8
4.2	Density of solids	8
4.2.1	Principle of density determination of solids	8
4.2.2	Determining the density of solids	9
5	Technical Data	12
5.1	Wires	12
5.2	Density of distilled water	12
6	Spare Parts and Accessories	13

1 Introduction

Thank you for choosing a METTLER TOLEDO density kit.

The density kit allows you to determine the density of solids. The kit is compatible with a wide range of advanced and standard balances. For a complete list of compatible balances, see [Compatibility table ▶ Page 6].

1.1 Further documents and information

▶ www.mt.com/lab-accessories

This document is available in other languages online.

▶ www.mt.com/Density-Kit-adv-std-RM

Search for documents:

▶ www.mt.com/library

For further questions, please contact your authorized METTLER TOLEDO dealer or service representative.

▶ www.mt.com/contact

1.2 Explanation of conventions and symbols used

Conventions and symbols

 **Note** For useful information about the product.



Refers to an external document.

Elements of instructions

In this manual, step-by-step instructions are presented as follows. The action steps are numbered and can contain prerequisites, intermediate results and results, as shown in the example. Sequences with less than two steps are not numbered.

■ Prerequisites that must be fulfilled before the individual steps can be executed.

1 Step 1

⇒ Intermediate result

2 Step 2

⇒ Result

1.3 Acronyms and abbreviations

Original term	Explanation
RM	Reference Manual

2 Safety information

A document named "Reference Manual" is available for this instrument.

- This Reference Manual contains a full description of the instrument and its use.
- Keep the Reference Manual for future reference.
- Include the Reference Manual if you transfer the instrument to other parties.

Only use the instrument according to the Reference Manual. If you do not use the instrument according to the Reference Manual or if it is modified, the safety of the instrument may be impaired and Mettler-Toledo GmbH assumes no liability.

2.1 Definitions of signal words and warning symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

Signal words

DANGER	A hazardous situation with high risk, resulting in death or severe injury if not avoided.
WARNING	A hazardous situation with medium risk, possibly resulting in death or severe injury if not avoided.
CAUTION	A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.
NOTICE	A hazardous situation with low risk, resulting in damage to the instrument, other material damage, malfunctions and erroneous results, or loss of data.

Warning symbols



Notice

2.2 Product-specific safety notes

Intended use

This instrument is intended to be used by trained staff. The density kit is designed to be installed on a range of METTLER TOLEDO analytical balances to determine the density of solids and liquids.

Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties.

Mettler-Toledo GmbH assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. Mettler-Toledo GmbH assumes that the instrument owner provides the necessary protective gear.

Safety notes




NOTICE

Damage to the instrument or malfunction due to the use of unsuitable parts

- Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

3 Installation and Putting into Operation




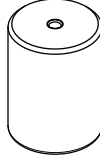


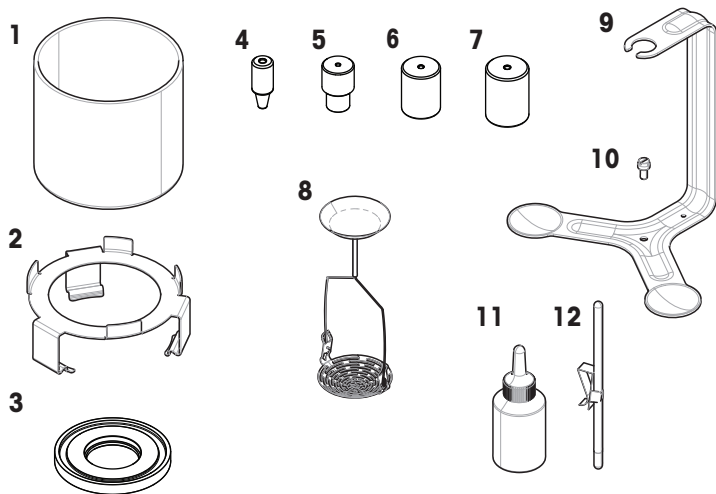
3.1 Scope of delivery



1	Beaker, \varnothing 85 mm, H 75 mm	8	Holder for sinking and floating solids
2	Beaker holder	9	Bracket
3	Adapter ring	10	Screw
4	Cone adapter	11	Wetting agent
5	Cone adapter	12	Thermometer, non-calibrated, with holder
6	Cone adapter (aluminum)		Reference Manual
7	Cone adapter (copper)		


3.2 Compatibility table

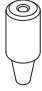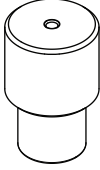Find the model, readability, and capacity of your balance on the type label.

Balance family	Readability	Cone adapter	Cone adapter	Cone adapter (aluminum)	Cone adapter (copper)	Adapter ring (depicted side facing up)
						
MX	0.01 mg	✓	–	–	–	
MX	0.1 mg	–	✓	–	–	
MX	1 mg	–	–	–	✓	
MR, MA	0.1 mg	–	✓	–	–	–
MA	0.01 mg	–	✓	–	–	
MR, MA	1 mg	–	–	✓	–	

3.3 Installing the density kit

- The proper compensation weight has been determined for the balance type, see [Compatibility table ▶ Page 6].
- 1 Disconnect the balance from the power supply.
- 2 Open the draft shield.
- 3 Remove the weighing pan.
- 4 Install the adapter ring (1) if needed. For correct orientation, see [Compatibility table ▶ Page 6].

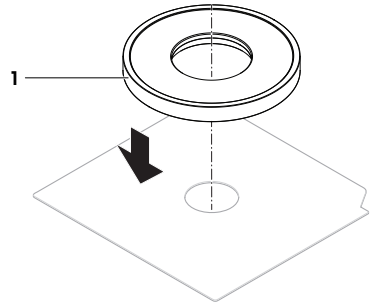


- 5 Assemble the bracket (2) and the appropriate compensation weight (3) with the provided screw (4).


Note

Depending on the balance model, the compensation weight in the illustration may look different.

- 6 Install the compensation weight and bracket on the weighing cell.

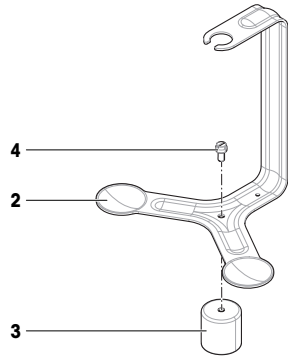

- 7 Place the beaker holder (5) on the balance (6). Make sure that the widest foot faces the front of the balance; the beaker holder must not be in contact with the bracket.
- 8 Reconnect the balance to the power supply.
- ➔ The density kit is ready to be used.

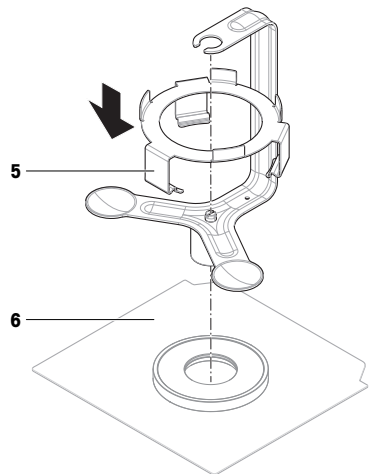

4 Operation

This section provides theoretical and practical information about how to determine the density of solids and liquids using the density kit. It also includes recommendations to improve the results' accuracy and precision. Depending on your balance model and software version, the main operation steps and necessary calculations might be implemented in the balance software; follow the instructions from the wizard for support.



For further information, consult the Reference Manual (RM) of your balance.

4.1 Principle of density determination

The density ρ is the quotient of the mass m and the volume V .

Density:

The international system of units specifies kg/m^3 as the unit of density. However, the unit g/cm^3 is better suited to laboratory purposes.

$$\rho = \frac{m}{V}$$

Density determinations are frequently performed by Archimedes' principle, which is also used with this density kit. **Archimedes' principle states that a body immersed partially or fully in fluid experiences a buoyant force acting upwards on it. The magnitude of this force is equivalent to the weight of the fluid displaced by the body.**

The procedure for the density determination by Archimedes' principle depends on whether the density of solids or liquids has to be determined.

4.2 Density of solids

4.2.1 Principle of density determination of solids

The density of a solid is determined with the aid of a liquid whose density ρ_0 is known. Water or ethanol are usually used as auxiliary liquids. The solid is weighed in air (A) and then in the auxiliary liquid (B). The density ρ can be calculated from the two weighing results.

Density:

$$\rho = \frac{A}{A-B} (\rho_0 - \rho_L) + \rho_L$$

Volume:

$$V = \alpha \frac{A-B}{\rho_0 - \rho_L}$$

Legend

A	Weight of sample in air
B	Weight of sample in the auxiliary liquid
ρ	Density of sample
ρ_0	Density of auxiliary liquid
ρ_L	Density of air (0.0012 g/cm^3)
V	Volume of sample
α	Balance correction factor (0.99985). It takes air buoyancy of the adjustment weight(s) into account.

4.2.2 Determining the density of solids



NOTICE

Erroneous results due to improper handling

The sample and the parts of the instrument immersed in liquid should be clean to obtain correct weighing results.


- Always use gloves when touching parts that are immersed in liquid.

To improve the accuracy of the results, take the following precautionary measures:



- The effect of temperature on liquids can cause density changes of order of magnitude 0.1 to 1 ‰ per °C. To obtain accurate results, take the temperature of the auxiliary liquid into account in all density determinations.
- Adhesion of the auxiliary liquid to the suspension wires of the basket causes an apparent weight increase of up to 3 mg. If the greatest possible accuracy is required, pour a few drops of the wetting agent in the auxiliary liquid.
- To avoid air bubbles:
 - Use enclosed wetting agent or organic liquids. The density change due to the addition of wetting agent can be neglected.
 - Remove grease traces from solvent-resistant solids.
 - Clean basket at regular intervals.
 - Remove adhering air bubbles with a fine brush.
 - Gently move the basket on first-time immersion to dislodge any air bubbles.

Preparing the density kit

- The draft shield doors are closed (if applicable).
- 1 Press →0← to zero the balance.
 - 2 Open the draft shield.
 - 3 Attach the supplied thermometer (1) to the beaker (2) and place the beaker on the pan (3).
 - 4 Fill the beaker with the auxiliary liquid. Add enough liquid to ensure that the solid is covered by at least 1 cm of liquid after immersion.

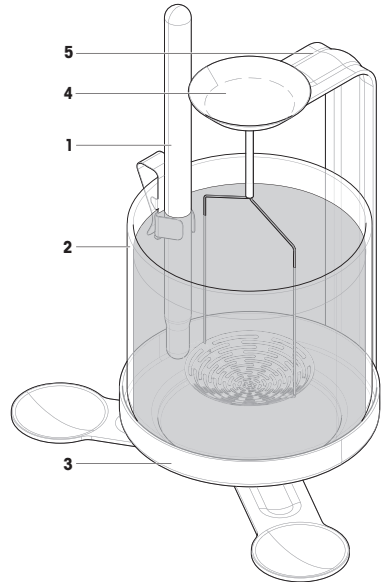


- 5 Carefully rotate the basket of the holder to the appropriate configuration: for sinking solids (4a) or floating (4b) solids.
- 6 Hang the holder (4) on the bracket (5). Make sure that no air bubbles adhere to the immersed part of the holder. Remove air bubbles either by moving the holder or by using a fine brush.
- 7 For floating solids: If the buoyancy of the solid is greater than the weight of the holder, the holder must be weighted by placing an additional weight (6) on the top weighing pan of the holder. This additional weight must remain on top of the weighing pan during the entire procedure.
- 8 Close the draft shield.
- 9 Press \rightarrow **T** \leftarrow to tare the balance.

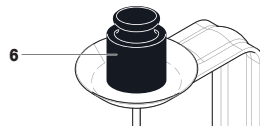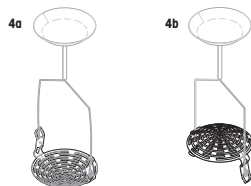



Weighing the sample in air

- 1 Open the draft shield.
- 2 Place the sample on one of the two weighing pans (1) of the bracket. If the weight of the sample is > 20 g, place it on the weighing pan at the top of the holder (2).
- 3 Wait until the balance has reached stability and note the displayed value (corresponding to "A" in the formula).
- 4 Open the draft shield.
- 5 Remove the sample from the weighing pan.

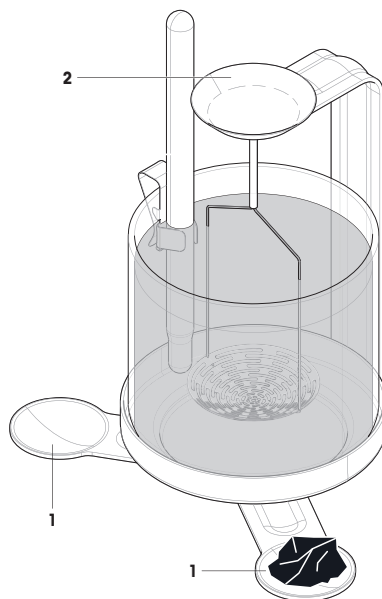

Weighing the sample in liquid

- 1 For sinking solids, place the sample in the holder's basket (A). For floating solids, place the sample under the holder's basket (B).
- 2 Ensure that no air bubbles adhere to the sample. Remove any air bubbles with a fine brush.
- 3 Close the draft shield.
- 4 Wait until the balance has reached stability and note the displayed value (corresponding to "B" in the formula).
- 5 Determine the density ρ of the solid according to the corresponding formula.

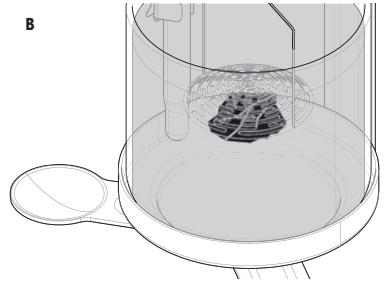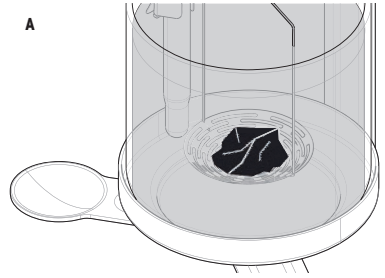

See also

- 🔗 Principle of density determination of solids ▶ Page 8

5 Technical Data

5.1 Wires


	Wire of holder
Material	Stainless steel
Diameter	0.6 mm

5.2 Density of distilled water

T/ °C	0	0.1	0.2	0.3	0.4	0.5	0.6	0.7	0.8	0.9
10	0.99970	0.99969	0.99968	0.99967	0.99966	0.99965	0.99964	0.99963	0.99962	0.99961
11	0.99960	0.99959	0.99958	0.99957	0.99956	0.99955	0.99954	0.99953	0.99952	0.99951
12	0.99950	0.99949	0.99947	0.99946	0.99945	0.99944	0.99943	0.99941	0.99940	0.99939
13	0.99938	0.99936	0.99935	0.99934	0.99933	0.99931	0.99930	0.99929	0.99927	0.99926
14	0.99924	0.99923	0.99922	0.99920	0.99919	0.99917	0.99916	0.99914	0.99913	0.99911
15	0.99910	0.99908	0.99907	0.99905	0.99904	0.99902	0.99901	0.99899	0.99897	0.99896
16	0.99894	0.99893	0.99891	0.99889	0.99888	0.99886	0.99884	0.99883	0.99881	0.99879
17	0.99877	0.99876	0.99874	0.99872	0.99870	0.99869	0.99867	0.99865	0.99863	0.99861
18	0.99859	0.99858	0.99856	0.99854	0.99852	0.99850	0.99848	0.99846	0.99844	0.99842
19	0.99840	0.99838	0.99836	0.99835	0.99833	0.99831	0.99828	0.99826	0.99824	0.99822
20	0.99820	0.99818	0.99816	0.99814	0.99812	0.99810	0.99808	0.99806	0.99803	0.99801
21	0.99799	0.99797	0.99795	0.99793	0.99790	0.99788	0.99786	0.99784	0.99781	0.99779
22	0.99777	0.99775	0.99772	0.99770	0.99768	0.99765	0.99763	0.99761	0.99758	0.99756
23	0.99754	0.99751	0.99749	0.99747	0.99744	0.99742	0.99739	0.99737	0.99734	0.99732
24	0.99730	0.99727	0.99725	0.99722	0.99720	0.99717	0.99715	0.99712	0.99709	0.99707
25	0.99704	0.99702	0.99699	0.99697	0.99694	0.99691	0.99689	0.99686	0.99684	0.99681
26	0.99678	0.99676	0.99673	0.99670	0.99667	0.99665	0.99662	0.99659	0.99657	0.99654
27	0.99651	0.99648	0.99646	0.99643	0.99640	0.99637	0.99634	0.99632	0.99629	0.99626
28	0.99623	0.99620	0.99617	0.99615	0.99612	0.99609	0.99606	0.99603	0.99600	0.99597
29	0.99594	0.99591	0.99588	0.99585	0.99582	0.99579	0.99577	0.99574	0.99571	0.99568
30	0.99564	0.99561	0.99558	0.99555	0.99552	0.99549	0.99546	0.99543	0.99540	0.99537

ISO 15212-1:1998 Oscillation-type density meters – Part 1: Laboratory instruments

6 Spare Parts and Accessories

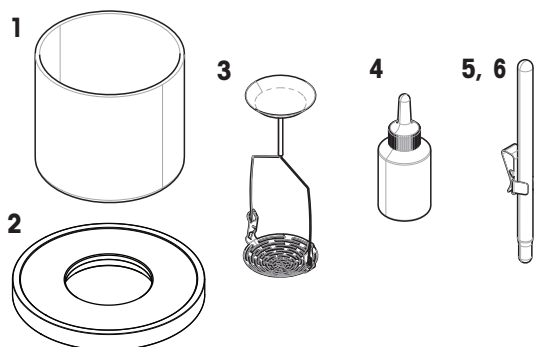

	Order no.	Designation	Remarks
1	30528711	Beaker, \varnothing 85 mm, H 75 mm	–
2	30881139	Adapter ring	–
3	30004746	Holder for sinking and floating solids	–
4	72409	Wetting agent	–
5	238767	Thermometer, non-calibrated	Including: holder
6	11132685	Thermometer, calibrated	Including: holder, calibration certificate

To protect your product's future:

METTLER TOLEDO Service assures the quality, measuring accuracy and preservation of value of this product for years to come.

Please request full details about our attractive terms of service.

► www.mt.com/service

www.mt.com/lab-accessories

For more information

Mettler-Toledo GmbH

Im Langacher 44
8606 Greifensee, Switzerland
www.mt.com/contact

Subject to technical changes.
© 10/2024 METTLER TOLEDO. All rights reserved.
30572182B en



30572182