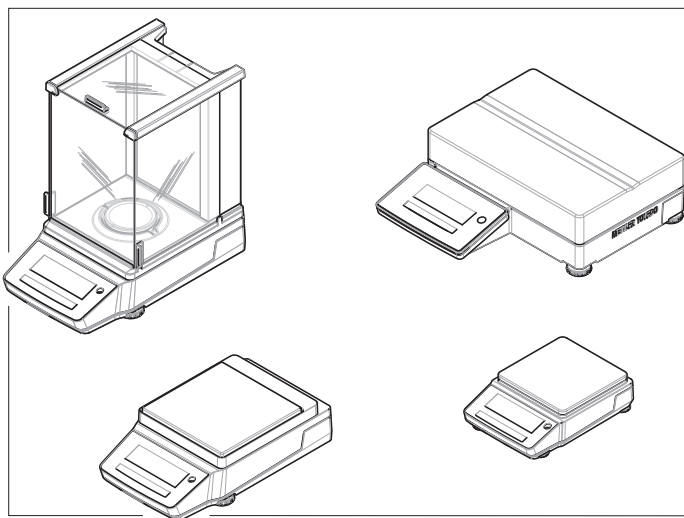
<b>English</b>	User Manual <b>Analytical and Precision Balances MA</b>
<b>Deutsch</b>	Benutzerhandbuch <b>Analysen- und Präzisionswaagen MA</b>
<b>Español</b>	Manual de usuario <b>Balanzas analíticas y de precisión MA</b>
<b>Français</b>	Guide de l'utilisateur <b>Balances d'analyse et de précision fine MA</b>
<b>简体中文</b>	用户手册 <b>分析与精密天平 MA</b>
<b>日本語</b>	ユーザマニュアル <b>分析天びん・上皿天びん MA</b>
<b>Čeština</b>	Návod k použití <b>Analytické a přesné váhy MA</b>
<b>Dansk</b>	Brugervejledning <b>Analyse- og præcisionsvægte MA</b>
<b>Hrvatski</b>	Korisnički priručnik <b>Analitičke i precizne vage MA</b>
<b>Magyar</b>	Felhasználói útmutató <b>Analitikai és precíziós mérlegek MA</b>

**METTLER TOLEDO**

► [www.mt.com/MA-RM](http://www.mt.com/MA-RM)

- en** For full information, always refer to the Reference Manual (RM).
- de** Ausführliche Informationen finden Sie stets im Referenzhandbuch (RM).
- es** Para obtener más información, consulte siempre el manual de referencia (RM).
- fr** Il convient de se reporter systématiquement au Manuel de référence (RM) pour obtenir des informations exhaustives.
- zh** 有关完整信息，请查阅《参考手册（RM）》。
- ja** すべての情報については、必ずリファレンスマニュアル（RM）を参照してください。
- cs** Podrobné informace naleznete vždy v referenční příručce (RM).
- da** Se altid referencemanualen (RM) for at få de komplette oplysninger.
- hr** Sve informacije uvijek možete pronaći u referentnom priručniku (RM).
- hu** A részletes információkat megtalálja a Referencia-kézikönyvben (RM).

---

User Manual **Analytical and Precision Balances**

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Magyar

# Table of Contents

<b>1</b>	<b>Introduction</b>	<b>3</b>
1.1	Document purpose .....	3
1.2	Further documents and information .....	3
1.3	Acronyms and abbreviations .....	3
<b>2</b>	<b>Safety Information</b>	<b>4</b>
2.1	Definition of signal words and warning symbols .....	4
2.2	Product-specific safety notes .....	4
<b>3</b>	<b>Design and Function</b>	<b>5</b>
3.1	Overview analytical balances .....	6
3.2	Overview precision balances, small .....	6
3.3	Overview precision balances, large .....	7
3.4	Overview precision balances, compact .....	7
3.5	Overview terminal .....	8
3.6	Overview interface connections .....	9
3.7	Components description .....	10
3.7.1	Draft shield .....	10
3.7.2	Weighing pan .....	10
3.7.3	Drip tray .....	10
3.7.4	Door handle .....	10
3.7.5	Leveling feet .....	11
3.7.6	Terminal .....	11
3.7.7	QuickLock for doors and front panel .....	11
3.7.8	Release button for back panel .....	11
3.8	User interface .....	12
3.8.1	Main sections at a glance .....	12
3.8.2	Main weighing screen .....	12
<b>4</b>	<b>Installation and Putting into Operation</b>	<b>15</b>
4.1	Selecting the location .....	15
4.2	Unpacking the balance .....	15
4.3	Installation .....	16
4.3.1	Balances with draft shield .....	16
4.3.2	Balances without draft shield .....	18
4.3.3	Balances, large .....	19
4.3.4	Balances, compact .....	19
4.4	Putting into operation .....	19
4.4.1	Connecting the balance .....	19
4.4.2	Switching on the balance .....	20
4.4.3	Leveling the balance .....	20
4.4.3.1	Leveling large balances .....	20
4.4.3.2	Leveling compact balances .....	21
4.4.3.3	Leveling all other balances .....	22
4.4.4	Performing an internal adjustment .....	22
4.4.5	Entering / Exiting standby mode .....	23
4.4.6	Entering / Exiting hibernation mode .....	23
4.4.7	Switching off the balance .....	23
4.5	Performing a simple weighing .....	23
4.5.1	Opening and closing draft shield doors .....	23
4.5.2	Zeroing the balance .....	23
4.5.3	Taring the balance .....	24
4.5.4	Performing a weighing .....	24

4.6	Using batteries (compact balance).....	24
4.6.1	Inserting or replacing batteries.....	24
4.7	Transporting, packing, and storing.....	25
4.7.1	Transporting the balance over short distances.....	25
4.7.2	Transporting the balance over long distances.....	25
4.7.3	Packing and storing.....	25
<hr/>		
<b>5</b>	<b>Maintenance</b>	<b>26</b>
5.1	Maintenance tasks.....	26
5.2	Cleaning.....	26
5.2.1	Disassembling for cleaning.....	26
5.2.1.1	Balances with draft shield.....	27
5.2.1.2	Balances without draft shield.....	29
5.2.1.3	Balances, large.....	29
5.2.1.4	Balances, compact.....	30
5.2.2	Cleaning the balance.....	30
5.2.3	Putting into operation after cleaning.....	31
5.3	Service.....	31
<hr/>		
<b>6</b>	<b>Technical Data</b>	<b>31</b>
6.1	General data.....	31
<hr/>		
<b>7</b>	<b>Disposal</b>	<b>32</b>
<hr/>		
<b>8</b>	<b>Compliance Information</b>	<b>32</b>

## 1 Introduction

Thank you for choosing a METTLER TOLEDO balance. The balance combines high performance with ease of use.

### EULA

The software in this product is licensed under the METTLER TOLEDO End User License Agreement (EULA) for Software.

When using this product you agree to the terms of the EULA.

▶ [www.mt.com/EULA](http://www.mt.com/EULA)

### 1.1 Document purpose

This User Manual provides brief instructions about the first steps to take with the instrument. This ensures safe and efficient handling. Personnel must have carefully read and understood this manual before performing any task.

### 1.2 Further documents and information

This document is available in other languages online.

▶ [www.mt.com/MA-UM](http://www.mt.com/MA-UM)

Product page:

▶ [www.mt.com/MA-balances](http://www.mt.com/MA-balances)

Instructions for cleaning a balance, "8 Steps to a Clean Balance":

▶ [www.mt.com/lab-cleaning-guide](http://www.mt.com/lab-cleaning-guide)

Search for software:

▶ [www.mt.com/labweighing-software-download](http://www.mt.com/labweighing-software-download)

Search for documents:

▶ [www.mt.com/library](http://www.mt.com/library)

For further questions, please contact your authorized METTLER TOLEDO dealer or service representative.

▶ [www.mt.com/contact](http://www.mt.com/contact)

### 1.3 Acronyms and abbreviations

Original term	Explanation
ABS/PC	Acrylonitrile butadiene styrene / polycarbonate (polymer blend)
AC	Alternating Current
ASTM	American Society for Testing and Materials
DC	Direct Current
EMC	Electromagnetic Compatibility
FCC	Federal Communications Commission
GWP	Good Weighing Practice
ID	Identification
IP	Ingress Protection
LPS	Limited Power Source
MT-SICS	METTLER TOLEDO Standard Interface Command Set
NA	Not Applicable
OIML	Organisation Internationale de Métrologie Légale

PA 12	Polyamid 12
PBT	Polybutylene terephthalate
PET	Polyethylene terephthalate
POM	Polyoxymethylene
RM	Reference Manual
SOP	Standard Operating Procedure
TDNR	Type Definition Number
TPE	Thermoplastic elastomer
UM	User Manual
USB	Universal Serial Bus
USP	United States Pharmacopeia

## 2 Safety Information

Two documents named "User Manual" and "Reference Manual" are available for this instrument.

- The User Manual is available online in various languages.
- A printed version of the User Manual is delivered with the instrument.
- The Reference Manual is available online. This manual contains a full description of the instrument and its use.
- Keep both documents for future reference.
- Include both documents if you transfer the instrument to other parties.

Only use the instrument according to the User Manual and the Reference Manual. If you do not use the instrument according to these documents or if the instrument is modified, the safety of the instrument may be impaired and Mettler-Toledo GmbH assumes no liability.

### 2.1 Definition of signal words and warning symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

#### Signal words

<b>DANGER</b>	A hazardous situation with high risk, resulting in death or severe injury if not avoided.
<b>WARNING</b>	A hazardous situation with medium risk, possibly resulting in death or severe injury if not avoided.
<b>CAUTION</b>	A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.
<b>NOTICE</b>	A hazardous situation with low risk, resulting in damage to the instrument, other material damage, malfunctions and erroneous results, or loss of data.

#### Warning symbols

General hazard

Notice

### 2.2 Product-specific safety notes

#### Intended use

This instrument is designed to be used by trained staff. The instrument is intended for weighing purposes.

Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

## Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties.

Mettler-Toledo GmbH assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. Mettler-Toledo GmbH assumes that the instrument owner provides the necessary protective gear.

## Safety notes

### **WARNING**

#### **Death or serious injury due to electric shock**

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.

### **NOTICE**

#### **Damage to the instrument or malfunction due to the use of unsuitable parts**

- Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

A list of spare parts and accessories can be found in the Reference Manual.

## 3 Design and Function

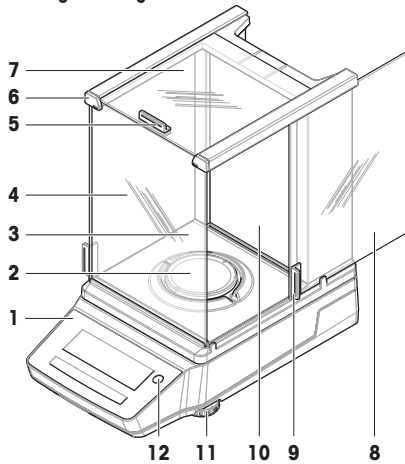
For further information, consult the Reference Manual (RM).

► [www.mt.com/MA-RM](http://www.mt.com/MA-RM)

### 3.1 Overview analytical balances


0.01 mg / 0.1 mg


1	Terminal	7	Top door, draft shield
2	Weighing pan	8	Side door, draft shield (right/left)
3	Drip tray	9	Handle, side door
4	Front panel, draft shield	10	Back panel, draft shield
5	Handle, top door	11	Leveling feet
6	QuickLock, front panel	12	Level indicator

### 3.2 Overview precision balances, small

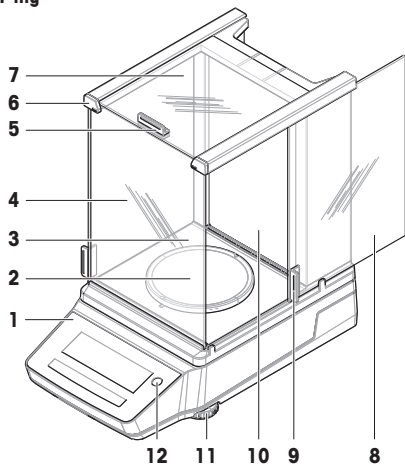1 mg

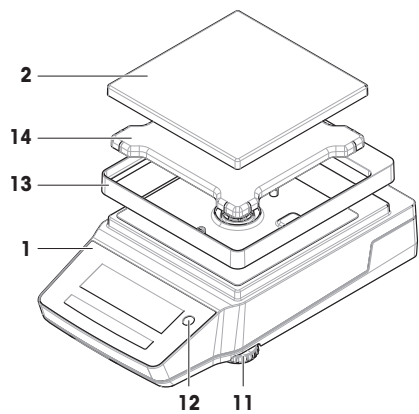
0.01 g / 0.1 g

1	Terminal	8	Side door, draft shield (right/left)
2	Weighing pan	9	Handle, side door
3	Drip tray	10	Back panel, draft shield
4	Front panel, draft shield	11	Leveling feet
5	Handle, top door	12	Level indicator
6	QuickLock, front panel	13	Draft-protection element
7	Top door, draft shield	14	Weighing pan support

### 3.3 Overview precision balances, large


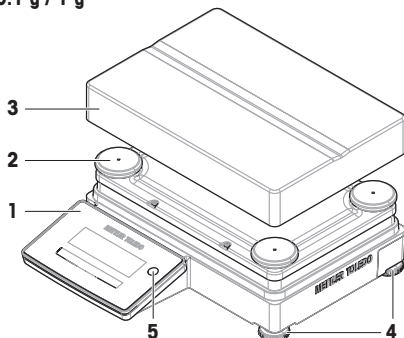
0.1 g / 1 g

1	Terminal	4	Leveling feet
2	Weighing pan support cap	5	Level indicator
3	Weighing pan		


### 3.4 Overview precision balances, compact

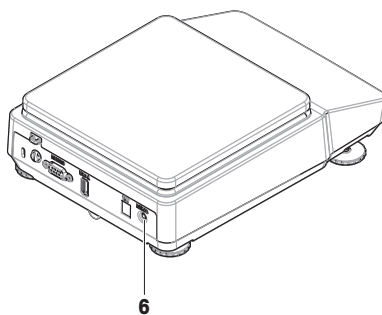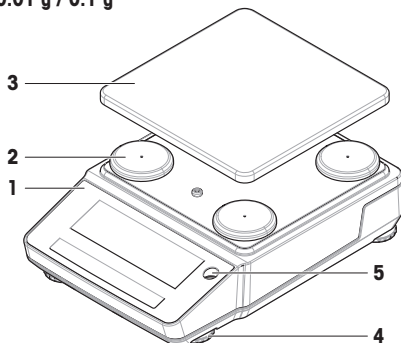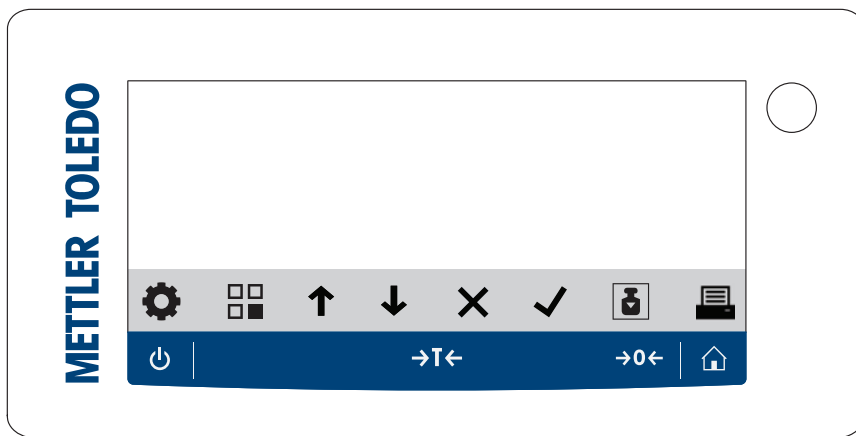0.01 g / 0.1 g

1	Terminal	4	Leveling feet
2	Weighing pan support cap	5	Level indicator
3	Weighing pan	6	Wake-up switch (battery mode)

### 3.5 Overview terminal

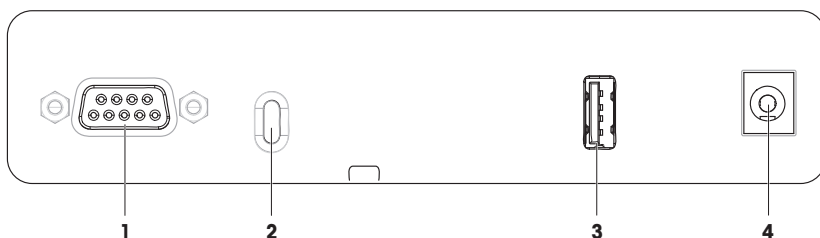
#### Terminal keys

Name	Description
 Standby	By tapping  , the balance is not completely switched off but goes into standby mode. To switch the balance completely off, it must be unplugged from the power supply. <b>Note</b> Do not disconnect the balance from the power supply unless the balance is not used for an extended period of time. After switching on the instrument, it must warm up before giving accurate results. <b>Note</b> Compact balances in battery mode only: By tapping  , the balance is completely switched off. There is no standby mode.
 Tare	Tares the balance. This function is used when the weighing process involves containers. After taring the balance, the screen shows <i>Net</i> which indicates that all displayed weights are net.
 Zero	Zeroes the balance. The balance must always be zeroed before starting the weighing process. After zeroing, the balance sets a new zero point.
 Home	To return from any settings menu level to the main weighing screen of the currently set or last used application.

## Operation keys

	Name	Description
	Settings	Accesses the settings menu.
	Applications	Accesses the applications menu.
	Previous / Up / Increase	<ul style="list-style-type: none"> <li>• Scrolls up in a list of topics.</li> <li>• Changes between unit 1, unit 2, and the application unit (if available).</li> <li>• Increases numbers.</li> <li>• Adds a captured weight within an application.</li> </ul>
	Next / Down / Decrease	<ul style="list-style-type: none"> <li>• Scrolls down in a list of topics.</li> <li>• Changes between unit 1, unit 2, and the application unit (if available).</li> <li>• Decreases numbers.</li> </ul>
	Cancel	<ul style="list-style-type: none"> <li>• Cancels a task or leaves the menu without saving.</li> <li>• Discards a sample in a workflow application.</li> </ul>
	Accept	<ul style="list-style-type: none"> <li>• Accesses the selected menu.</li> <li>• Starts the selected application.</li> <li>• Confirms the entry.</li> </ul>
	Adjustment	Executes the predefined adjustment procedure.
	Print / Transfer	<ul style="list-style-type: none"> <li>• Prints the displayed value.</li> <li>• Transfers data to the predefined target location.</li> </ul>

### 3.6 Overview interface connections


<b>1</b>	RS232C serial interface	<b>3</b>	USB-A port
<b>2</b>	Slot for anti-theft cable	<b>4</b>	Socket for AC/DC adapter

### 3.7 Components description

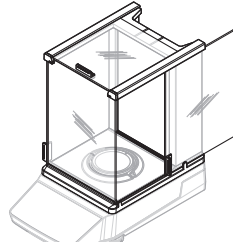
#### 3.7.1 Draft shield

The draft shield protects the weighing area against environmental impacts like drafts or moisture. The side doors and the top door can be opened manually.


#### 3.7.2 Weighing pan

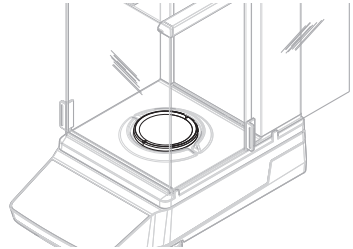The weighing pan is the load receptor that serves to accommodate the weighing item.


#### 3.7.3 Drip tray

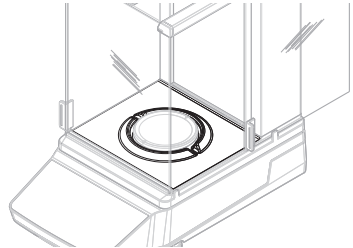The drip tray is positioned below the weighing pan. The primary purpose of the drip tray is to ensure quick cleaning of the balance.

#### 3.7.4 Door handle

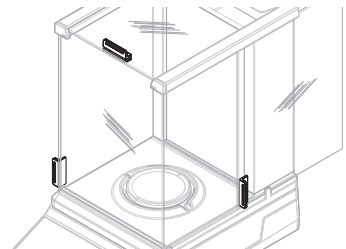The door handles are mounted on the draft shield doors. The handles are used to manually open the side doors and the top door of the draft shield.


### 3.7.5 Leveling feet

The balance stands on height-adjustable feet. These feet are used to level the balance.


**Note**

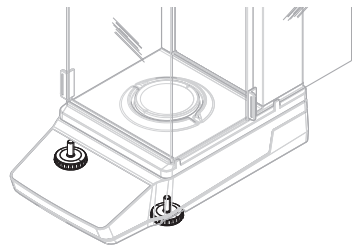Large balances and compact balances have four leveling feet.


### 3.7.6 Terminal

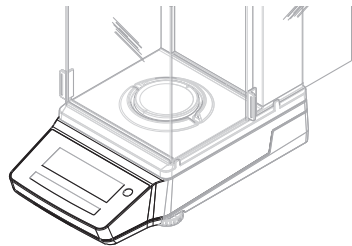The terminal is integrated into the balance and has an LCD display. The terminal and the platform are protected by a replaceable cover.


### 3.7.7 QuickLock for doors and front panel

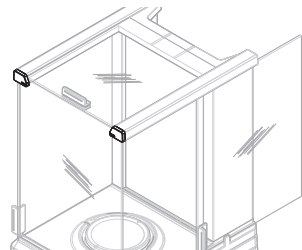Depending on the position, the QuickLock is used to lock/unlock the top door, the side doors, and the front panel of the draft shield.

### 3.7.8 Release button for back panel

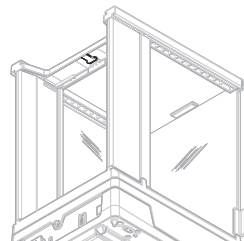The release button is used to lock/unlock the back panel of the draft shield.

### 3.8 User interface

#### 3.8.1 Main sections at a glance

The main weighing screen (1) displays weighing results and provides feedback on the status of the balance. The operation keys (2) are the main navigation element where all the menus and settings can be accessed and selected.

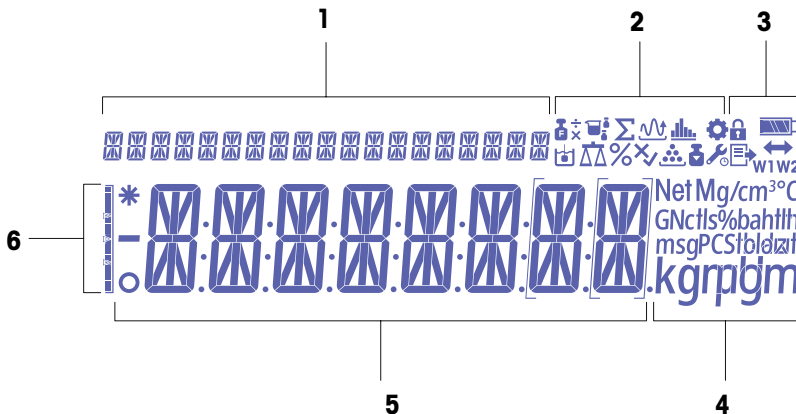
	Name	Description
1	Main weighing screen	Displays the weighing results and provides information on the status of the balance.

#### See also

[Main weighing screen](#) ▶ Page 12

#### 3.8.2 Main weighing screen

1	Coach text	4	Unit icons
2	Application icons	5	Weight value / menu label / parameter label / parameter value
3	Status icons	6	Weighing aid

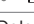
#### Application icons

	Name	Description
	Application "Weighing"	The application "Weighing" is used for simple weighing tasks.
	Application "Piece counting"	The application "Piece counting" allows you to determine the number of pieces put on the weighing pan.

	Name	Description
	Application "Percent weighing"	"Percent weighing" allows a sample weight to be checked as a percentage of a reference target weight
	Application "Formulation"	The application "Formulation" (Net total) has several purposes. <ul style="list-style-type: none"> <li>• Weigh in (add and store) up to 999 individual component weights and display the total. If a printer is connected, the component weights are printed individually and as a total.</li> <li>• Tare and store up to 999 container weights and display the total. If a printer is connected, the tare weights are printed individually and as a total.</li> <li>• Fill up the sum of all component net weight values by adding a further component to a higher value.</li> </ul>
	Application "Dynamic weighing"	The application "Dynamic weighing" allows you to determine the weights of unstable samples or to determine weights under unstable ambient conditions. The balance calculates the weight as the average of a number of weighing operations over a defined time.
	Application "Density"	The application "Density" allows you to determine the density of solid bodies. Determination of the density uses "Archimedes' principle" according to which a body immersed in a fluid undergoes an apparent loss in weight which is equal to the weight of the fluid it displaces.
	Application "Check weighing"	The application "Check weighing" allows you to check the deviation of a sample weight within a tolerance limit to a reference target weight.
	Application "Factor weighing"	The application "Factor weighing" multiplies a predefined factor by the weight value (in grams) or divides the weight value by a predefined factor. Depending on the method used, one the following equation applies: <ul style="list-style-type: none"> <li>• result = weight / factor</li> <li>• result = weight × factor</li> <li>• result = weight + factor</li> <li>• result = weight – factor</li> </ul> The result is rounded to a predefined number of decimal places.
	Application "Statistics"	The application "Statistics" allows you to generate statistics of a series of weighing values. 1 to 999 values are possible.
	Application "Totaling"	The application "Totaling" allows you to weigh in different samples to add their weight values and to totalize them. 1 to 999 samples are possible.

### Status icons

	Name	Description
	In the menu "Settings"	The Menu settings is open and an entry of this menu can be selected.
	Access protection on	Menu settings are locked and cannot be adapted.
	Adjustments started	The internal or external adjustment of the balance has started.
	Service reminder	<ul style="list-style-type: none"> <li>• Icon is blinking: next service is due.</li> <li>• Icon lights up constantly: service mode is active.</li> </ul>
	Publishing	Data is being published. If this icon is displayed, it is not possible to start a new data transfer.

	Name	Description
	Status of battery charge	Compact balances only: Indicates the level of the battery charge. <ul style="list-style-type: none"> <li>: Battery full</li> <li>: Battery 3/4 full</li> <li>: Battery 1/2 full</li> <li>: Battery 1/4 full</li> <li>: Battery discharged</li> </ul>
	Connectivity	Data connection to a USB device.
<b>W1</b>	Weighing range 1	Weighing range 1 is active.
<b>W2</b>	Weighing range 2	Weighing range 2 is active.
<b>Net</b>	Net weight values	"Net" indicates that all weight values displayed are net values.
<b>G</b>	Gross weight values	Gross weight values are displayed.
<b>M</b>	Stored value (Memory)	A stored value from the memory of the balance is displayed.
<b>X</b>	Failed operation	An operation has not been completed successfully.

#### Label Icons

	Name	Description
<b>—</b>	Negative value	Indicates that the displayed values are negative.
<b>*</b>	Calculated value	Indicates that the displayed value is calculated.
<b>o</b>	Unstable value	Indicates the displayed value is unstable, meaning that it changes over time.
<b>[ ]</b>	Uncertified digits	Brackets indicate uncertified digits (approved balance models only). <ul style="list-style-type: none"> <li><b>[ ]</b>: First decimal place.</li> <li><b>[ ]</b>: First decimal place for dual range balance.</li> </ul>

## 4 Installation and Putting into Operation

### 4.1 Selecting the location

A balance is a sensitive precision instrument. The location where it is placed will have a profound effect on the accuracy of the weighing results.

#### Requirements of the location

Place indoors on stable table

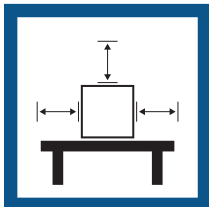
Ensure sufficient spacing

Level the instrument

Provide adequate lighting

Avoid direct sunlight



Avoid vibrations

Avoid strong drafts

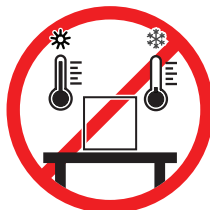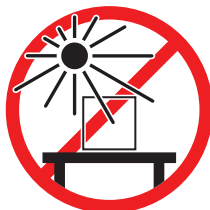Avoid temperature fluctuations

Take into account the environmental conditions. See "Technical Data".

Sufficient spacing for balances: > 15 cm all around the instrument


### 4.2 Unpacking the balance

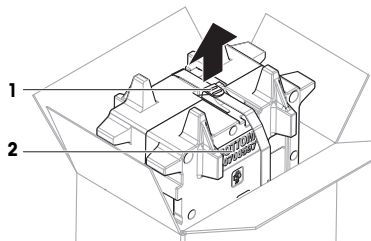Check the package, the packaging elements and the delivered components for damages. If any components are damaged, please contact your METTLER TOLEDO service representative.


#### **Note**

Depending on the balance model, the packaging elements and the components may look different.


- 1 Open the box and lift the package out using the lifting strap (1).
- 2 Place the package on a level surface with the inscription **BOTTOM (2)** facing downwards.

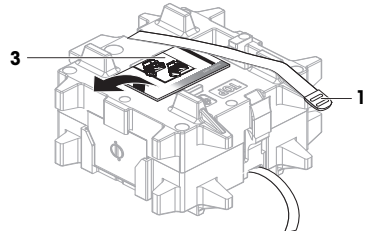
- 3 Open the lifting strap (1) and remove the User Manual (3).

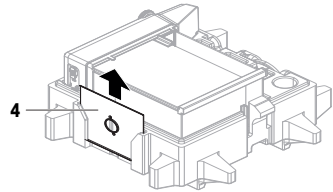- 4 Remove the upper part of the package and unpack the drip tray (4).

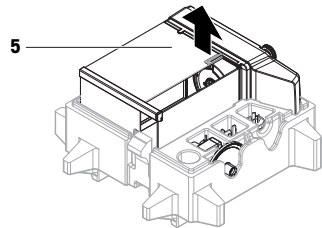- 5 Carefully unpack the balance (5) and all other items.
- 6 Remove the protective bag.
- 7 Keep the protective cover installed on the platform and on the terminal.
- 8 Store all parts of the packaging in a safe place for future use.
  - ➔ The balance is ready for installation.

## 4.3 Installation

### Note

Depending on the balance model, the components may look different.

### 4.3.1 Balances with draft shield

### CAUTION

#### Injury due to sharp objects or broken glass

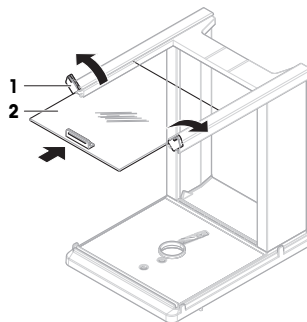
Instrument components, e.g., glass, can break and lead to injuries.

- Always proceed with focus and care.


**Note**

Skip steps 1 – 5 when unpacking the balance for the first time, or if the draft shield is already assembled and mounted on the platform.


- 1 Assemble the draft shield: Turn the QuickLock (1, right, left) and slide in the top door (2).

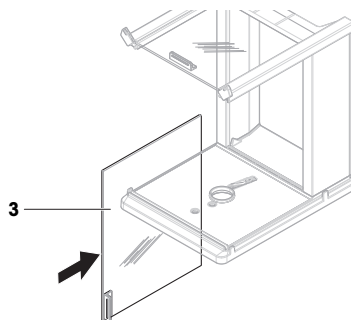
- 2 Slide in the side door (3) (right, left).

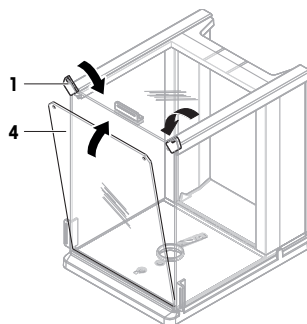
- 3 Attach the front panel (4), then turn the QuickLock (1, right, left) to hold the panel in place.

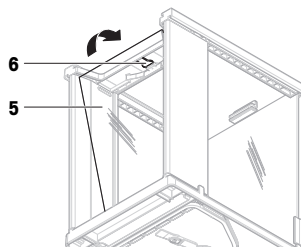- 4 Attach the back panel (5). Make sure the release button (6) snaps in.

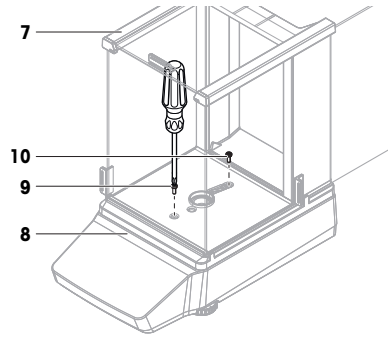
- Secure the draft shield (7) to the platform (8) by fixing the front screw (9) and the rear screw (10) with a Phillips screwdriver.


- Fully open the side door (11).
- Insert the drip tray (12).

**Note**

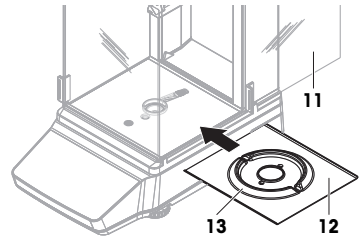The ring-shaped elevation (13) must point upwards. It serves as draft protection.

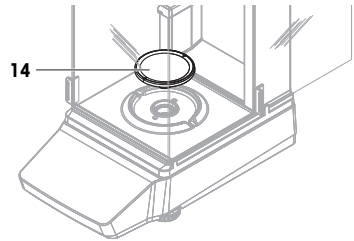
- Install weighing pan (14).
  - ➔ The balance is ready for use.

### 4.3.2 Balances without draft shield

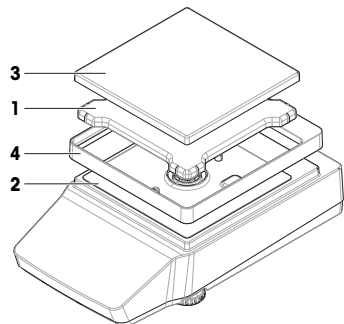- Place the weighing pan support (1) on top of the platform (2).
  - Note**  
To protect your balance, keep the protective cover installed on the platform (2).
- Place the weighing pan (3) on top of the weighing pan support (1).
- Place the draft-protection element (4) on top of the platform (2).
  - ➔ The balance is ready for use.


### 4.3.3 Balances, large

- 1 Remove the transport protection (1).
- 2 Place the weighing pan (2) on top of the support caps (3).

**Note**

To protect your balance, keep the protective cover installed on the terminal (4).

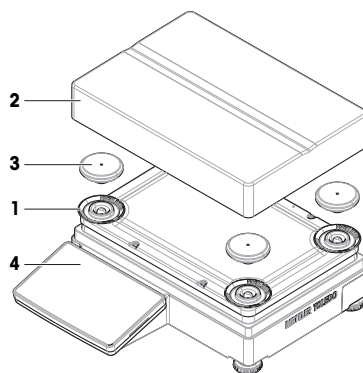⇒ The balance is ready for use.


### 4.3.4 Balances, compact

- Place the weighing pan (1) on top of the support caps (2).

**Note**

To protect your balance, keep the protective cover installed on the platform (3).

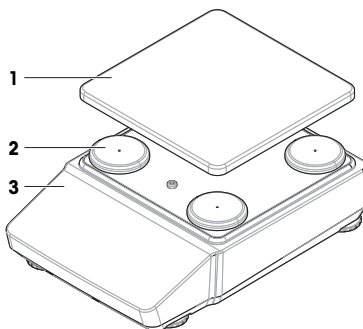⇒ The balance is ready for use.

## 4.4 Putting into operation

### 4.4.1 Connecting the balance

**WARNING**

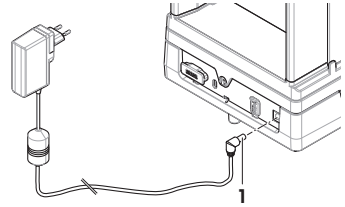
**Death or serious injury due to electric shock**

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.

- 1 Install the cables in such a way that they cannot be damaged or interfere with operation.

- 2 Insert the plug of the AC/DC adapter (1) into the power socket of the instrument.
- 3 Insert the plug of the power cable into a grounded power outlet that is easily accessible.
  - ➔ The balance automatically switches on.

**Note**

Do not connect the instrument to a power outlet controlled by a switch. After switching on the instrument, it must warm up before giving accurate results.

**See also**

🔗 General data ▶ Page 31

#### 4.4.2 Switching on the balance

When connected to the power supply, the balance automatically switches on.

##### Acclimatization and warm up

Before the balance gives reliable results, it must:

- acclimatize to the room temperature
- warm up by being connected to the power supply

The acclimatization time and warm-up time for balances are available in "General data".

**Note**

When the balance is exiting standby, it is ready immediately.

**See also**

🔗 General data ▶ Page 31

🔗 Entering / Exiting standby mode ▶ Page 23

🔗 Entering / Exiting hibernation mode ▶ Page 23

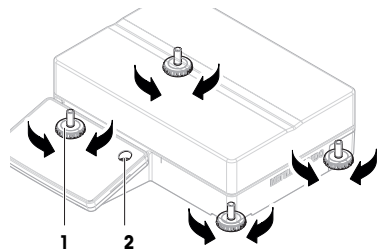
🔗 Switching off the balance ▶ Page 23

#### 4.4.3 Leveling the balance

Exact horizontal and stable positioning are essential for repeatable and accurate weighing results.

##### 4.4.3.1 Leveling large balances

- 1 Screw in all leveling feet fully.
- 2 Screw out all leveling feet two to three turns except for the rear left leveling foot.
- 3 Turn the leveling feet (1) until the air bubble is in the center of the level indicator (2). Proceed as shown in the following example.
- 4 Screw out the rear left leveling foot until it makes contact with the table surface.
  - ➔ The balance is leveled and supported by all four leveling feet.


**Example**

Air bubble at 12 o'clock:

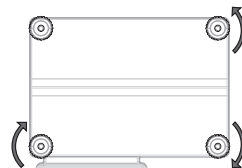
Turn both front feet clockwise.  
Turn rear right foot counterclockwise.

Air bubble at 3 o'clock:

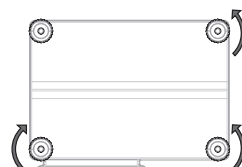
Turn left front foot clockwise, turn right front foot counterclockwise.  
Turn rear right foot counterclockwise.

Air bubble at 6 o'clock:

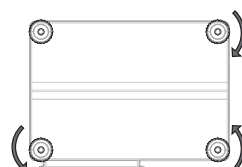
Turn both front feet counterclockwise.  
Turn rear right foot clockwise.


Air bubble at 9 o'clock:

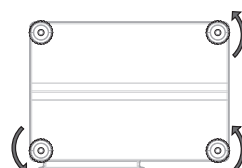Turn both front feet counterclockwise.  
Turn rear right foot counterclockwise.

**4.4.3.2 Leveling compact balances**

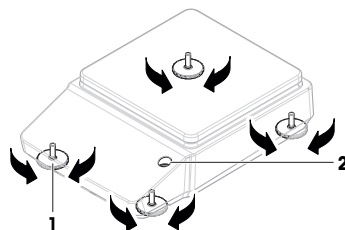- 1 Screw in both rear leveling feet fully.
- 2 Screw out both front leveling feet two to three turns.
- 3 Turn the front leveling feet (1) until the air bubble is in the center of the level indicator (2). Proceed as shown in the following example.
- 4 Screw out both rear leveling feet until they make contact with the table surface.
  - ➔ The balance is leveled and supported by all four leveling feet.

### Example

Air bubble at 12 o'clock:

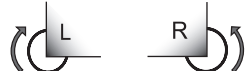
Turn both feet clockwise.

Air bubble at 3 o'clock:

Turn left foot clockwise, turn right foot counterclockwise.

Air bubble at 6 o'clock:

Turn both feet counterclockwise.

Air bubble at 9 o'clock:


Turn left foot counterclockwise, turn right foot clockwise.

#### 4.4.3.3 Leveling all other balances

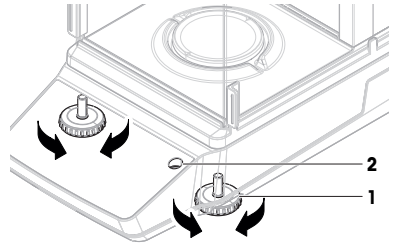- Turn the leveling feet (1) until the air bubble is in the center of the level indicator (2).


### Example

Air bubble at 12 o'clock:

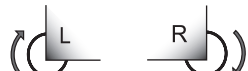
Turn both feet clockwise.

Air bubble at 3 o'clock:

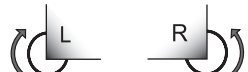Turn left foot clockwise, turn right foot counterclockwise.

Air bubble at 6 o'clock:

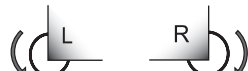
Turn both feet counterclockwise.

Air bubble at 9 o'clock:


Turn left foot counterclockwise, turn right foot clockwise.

#### 4.4.4 Performing an internal adjustment

- 1 On the main weighing screen, press .  
➔ ADJ.INT appears.
- 2 Press .
- ➔ The adjustment is executed.  
➔ The adjustment results appear.
- 3 Press .
- ➔ The balance is ready.
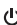
#### 4.4.5 Entering / Exiting standby mode

- 1 To enter standby mode, press  for less than 2 s.
  - ➔ The brightness of the display is reduced, information on the display is visible. The balance is still switched on.
- 2 To exit standby mode, press .
  - ➔ The display is turned on.

#### 4.4.6 Entering / Exiting hibernation mode

##### Note

Only relevant for compact balances in battery mode.

- 1 To enter hibernation mode, press  for less than 2 s.
  - ➔ The display is turned off. The balance is in hibernation mode.
- 2 To exit hibernation mode, press the wake-up switch on the rear side of the balance.
  - ➔ The display is turned on.

#### 4.4.7 Switching off the balance

To completely switch off the balance, it must be disconnected from the power supply. By pressing , the balance goes only into standby mode.

##### Note

When the balance has been completely switched off for some time, it must warm up before it can be used.

##### See also

-  Switching on the balance ▶ Page 20
-  Entering / Exiting standby mode ▶ Page 23
-  Entering / Exiting hibernation mode ▶ Page 23

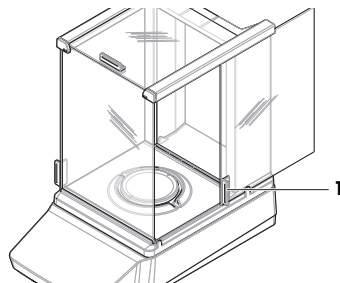
### 4.5 Performing a simple weighing

##### Note

A balance with draft shield is used to explain the procedure. For balances without a draft shield, skip the instructions steps concerning the draft shield.

#### 4.5.1 Opening and closing draft shield doors

- Open the door manually with the door handle (1).

#### 4.5.2 Zeroing the balance

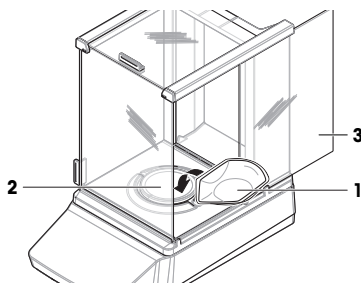
- 1 Open the draft shield.
- 2 Clear the weighing pan.
- 3 Close the draft shield.
- 4 Press  to zero the balance.
  - ➔ The balance is zeroed.

### 4.5.3 Taring the balance

If a sample vessel is used, the balance must be tared.

- The balance is zeroed.

  - 1 Place the sample vessel (1) on the weighing pan (2).
  - 2 Close the draft shield door (3).
  - 3 Press →T← to tare the balance.
    - ➔ The balance is tared. The icon  appears.

### 4.5.4 Performing a weighing

- 1 Open the draft shield.
- 2 Place the weighing object into the sample vessel.
- 3 Close the draft shield.
- 4 Wait until the instability detector  disappears.
  - ➔ The result is displayed.
- 5 Optional, if a printer is connected: Press  to print the weighing result.

## 4.6 Using batteries (compact balance)

The balance can also be operated with batteries. Under normal operating conditions, the balance runs independently of the AC power supply for about 8 to 15 hours (using alkaline batteries).

Immediately after an interruption of the AC power supply, the balance automatically switches to battery operation. This can happen, for example, by pulling the power plug, or in the event of a power failure. After the AC power supply is restored, the balance automatically switches back to AC operation.

The balance uses eight standard AA batteries (LR6). Alkaline batteries are preferred.

Rechargeable batteries can be used. Charging batteries inside the balance is **not** possible.

When the balance is operated with batteries, the battery symbol in the display lights up. The number of segments that are shown in the battery symbol indicate the charge level. When the batteries are almost completely discharged, the battery symbol flashes.

### See also

 Main weighing screen ▶ Page 12

### 4.6.1 Inserting or replacing batteries

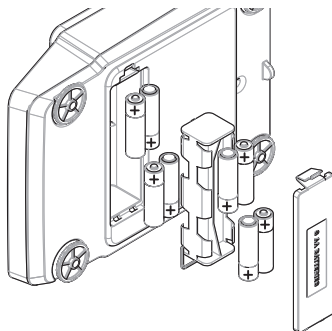
- Read and follow all warnings and instructions supplied by the battery manufacturer.
- Do not mix different types or brands of batteries. Performance of batteries varies depending on the manufacturer.
- Remove the batteries from the balance if the balance is not used for a long period of time.
- Batteries must be disposed of properly, according to local regulations.

Proceed as follows:

- The balance is switched off.
- The weighing pan is removed.

  - 1 Turn the balance carefully on its side.

- 2 Open and remove the battery-chamber cover.
- 3 Insert / replace the batteries with the correct polarity as shown in the battery holder.
- 4 Insert and close the battery-chamber cover.
- 5 Turn the balance carefully to its normal position.
- 6 Reinstall all components in the reverse order.


## 4.7 Transporting, packing, and storing

### 4.7.1 Transporting the balance over short distances

- 1 Disconnect the AC/DC adapter and unplug all interface cables.
- 2 Hold the balance with both hands and carry it in horizontal position to the target location. Consider the requirements of the location.

If you want to put the balance into operation, proceed as follows:

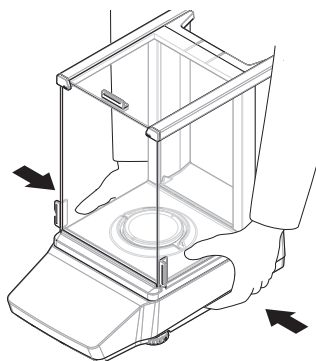- 1 Connect in reverse order.
- 2 Give the balance sufficient time to warm up.
- 3 Level the balance.
- 4 Perform an internal adjustment.

#### See also

- 🔗 Selecting the location ▶ Page 15
- 🔗 Switching on the balance ▶ Page 20
- 🔗 Leveling the balance ▶ Page 20
- 🔗 Performing an internal adjustment ▶ Page 22

### 4.7.2 Transporting the balance over long distances

METTLER TOLEDO recommends using the original packaging for transportation or shipment of the balance or balance components over long distances. The elements of the original packaging are developed specifically for the balance and its components and ensure maximum protection during transportation.

#### See also

- 🔗 Unpacking the balance ▶ Page 15

### 4.7.3 Packing and storing

#### Packing the balance

Store all parts of packaging in a safe place. The elements of the original packaging are developed specifically for the balance and its components, and ensures maximum protection during transportation and storage.

#### Storing the balance

Only store the balance under the following conditions:

- Indoor and in the original packaging

- According to the environmental conditions, see "Technical Data"

**Note**

When storing for longer than two weeks, the battery (capacitor) may become empty (only date and time get lost).

**See also**

 Technical Data ▶ Page 31

## 5 Maintenance

To guarantee the functionality of the balance and the accuracy of the weighing results, a number of maintenance actions must be performed by the user.

For further information, consult the Reference Manual (RM).

[www.mt.com/MA-RM](http://www.mt.com/MA-RM)

### 5.1 Maintenance tasks

Maintenance action	Recommended interval	Remarks
Performing an internal adjustment	<ul style="list-style-type: none"> <li>• Daily</li> <li>• After cleaning</li> <li>• After leveling</li> <li>• After changing the location</li> </ul>	see "Performing an internal adjustment"
Performing routine tests (eccentricity test, repeatability test, sensitivity test). METTLER TOLEDO recommends to at least perform a sensitivity test.	<ul style="list-style-type: none"> <li>• After cleaning</li> <li>• After assembling the balance</li> <li>• After a software update</li> <li>• Depending on your internal regulations (SOP)</li> </ul>	see "Tests" in the Reference Manual
Cleaning	<ul style="list-style-type: none"> <li>• After every use</li> <li>• Depending on the degree of pollution</li> <li>• Depending on your internal regulations (SOP)</li> </ul>	see "Cleaning"
Updating the software	<ul style="list-style-type: none"> <li>• Depending on your internal regulations (SOP).</li> <li>• After a new software release.</li> </ul>	see "Software update" in the Reference Manual

**See also**

 Performing an internal adjustment ▶ Page 22

 Cleaning ▶ Page 26

### 5.2 Cleaning

#### 5.2.1 Disassembling for cleaning

**Note**

Depending on the balance model, the components may look different.

**Note**

In most cases, it is not necessary to remove the protective cover to clean the balance.

### 5.2.1.1 Balances with draft shield

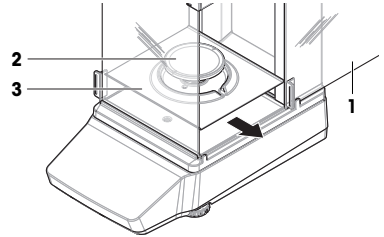
#### CAUTION


##### Injury due to sharp objects or broken glass

- Instrument components, e.g., glass, can break and lead to injuries.
- Always proceed with focus and care.


- 1 Fully open the side door (1).
- 2 Remove the weighing pan (2) and the drip tray (3).

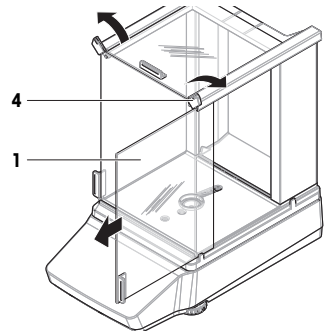
- 3 Turn the quick lock (4) and pull the side door (1) towards the front to remove it (right, left).

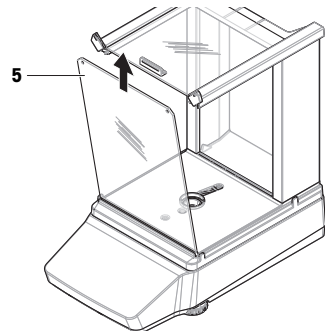- 4 Tilt the front panel (5) towards the front and lift it upwards to remove it.

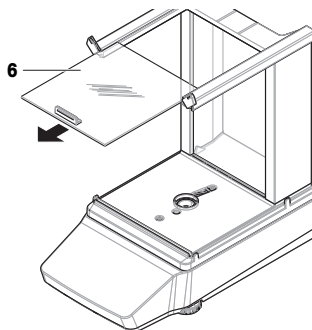
- 5 Pull the top door (6) towards the front to remove it.

- 6 Press the release button (7) and tilt the back panel (8) to remove it.


**Note**

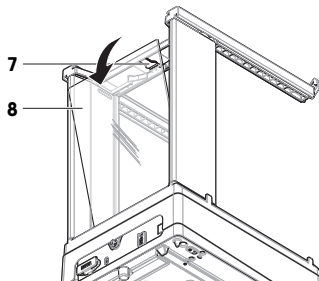Optional, if required: Remove the protective cover for cleaning as described below.


- 7 Remove the front screw (9) and the rear screw (10) with a Phillips screwdriver.

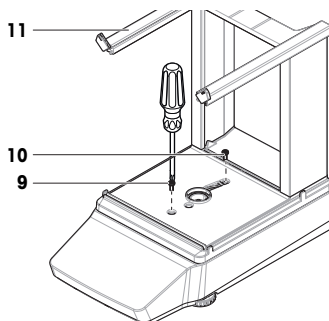- 8 Remove the draft shield (11).

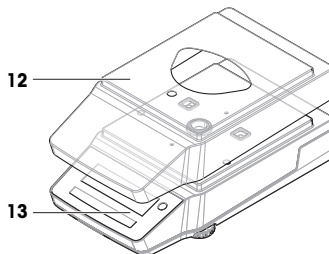- 9 Remove the protective cover (12) from the platform (13).


### 5.2.1.2 Balances without draft shield


- 1 Remove the draft-protection element (1).
- 2 Remove the weighing pan (2).
- 3 Remove the weighing pan support (3).

**Note**

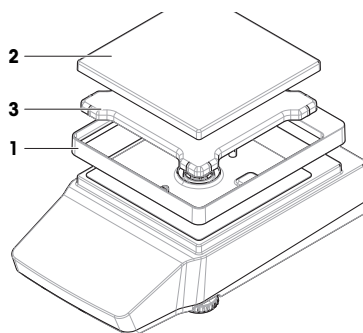Optional, if required: Remove the protective cover for cleaning as described below.

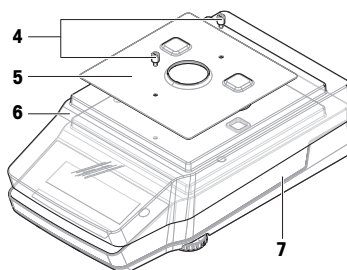
- 4 Remove the screws (4) to remove the EMC plate (5).
- 5 Remove the protective cover (6) from the platform (7).

### 5.2.1.3 Balances, large

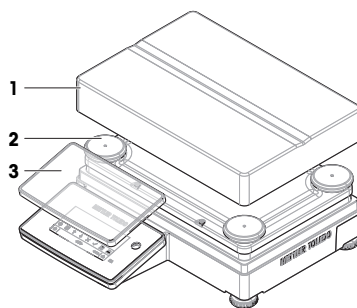- 1 Remove the weighing pan (1).
- 2 Remove the support caps (2).
- 3 Optional, if required: Remove the protective cover (3) for cleaning.


### 5.2.1.4 Balances, compact

- 1 Remove the weighing pan (1).
- 2 Remove the support caps (2).

**Note**

Optional, if required: Remove the protective cover for cleaning as described below.

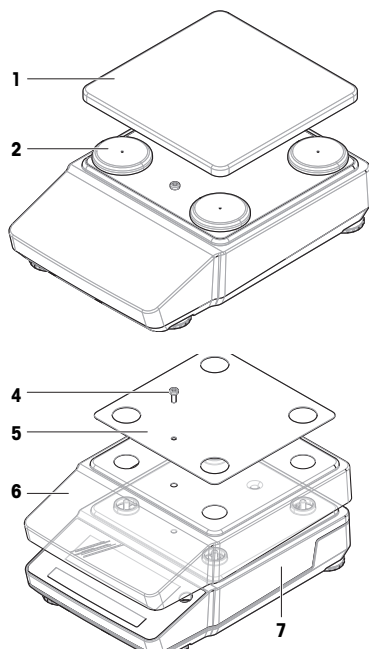- 3 Remove the screw (4) to remove the EMC plate (5).
- 4 Remove the protective cover (6) from the platform (7).

### 5.2.2 Cleaning the balance

#### NOTICE

##### Damage to the instrument due to inappropriate cleaning methods

If liquid enters the housing, it can damage the instrument. The surface of the instrument can be damaged by certain cleaning agents, solvents, or abrasives.

- 1 Do not spray or pour liquid on the instrument.
- 2 Only use the cleaning agents specified in the Reference Manual (RM) of the instrument or the guide "8 Steps to a Clean Balance".
- 3 Only use a lightly moistened, lint-free cloth or a tissue to clean the instrument.
- 4 Wipe off any spills immediately.

For further information on cleaning a balance, consult "8 Steps to a Clean Balance".

► [www.mt.com/lab-cleaning-guide](http://www.mt.com/lab-cleaning-guide)

#### Cleaning around the balance

- Remove any dirt or dust around the balance and avoid further contaminations.

#### Cleaning the terminal

- Clean the terminal with a damp cloth or a tissue and a mild cleaning agent.

### Cleaning the removable parts

- Clean the removed part with a damp cloth or a tissue and a mild cleaning agent or clean in a dishwasher up to 80 °C.

### Cleaning the balance

- 1 Disconnect the balance from the AC/DC adapter.
- 2 Use a lint-free cloth moistened with a mild cleaning agent to clean the surface of the balance.
- 3 Remove powder or dust with a disposable tissue first.
- 4 Remove sticky substances with a damp lint-free cloth and a mild solvent, e.g., isopropanol or ethanol 70%.

### 5.2.3 Putting into operation after cleaning

- 1 Reassemble the balance.
- 2 Check that the draft shield doors (top, sides) open and close normally (if applicable).
- 3 Reconnect the balance to the AC/DC adapter.
- 4 Check the level bubble, level the balance if necessary.
- 5 Respect the warm-up time specified in the "Technical Data".
- 6 Perform an adjustment.
- 7 Perform a routine test according to the internal regulations of your company.
- 8 Press **→0←** to zero the balance.
  - ⇒ The balance is ready for use.

#### See also

- 🔗 Leveling the balance ▶ Page 20
- 🔗 Technical Data ▶ Page 31
- 🔗 Performing an internal adjustment ▶ Page 22

## 5.3 Service

Regular servicing by an authorized service technician ensures reliability for years to come. Contact your METTLER TOLEDO representative for details about the available service options.

## 6 Technical Data

### 6.1 General data

#### Power supply

AC/DC adapter:	Input: 100 – 240 V AC ± 10%, 50 – 60 Hz, 0.5 A Output: 12 V DC, 1 A, LPS
Balance power consumption:	12 V DC, 0.5 A
Polarity:	

#### Protection and standards

Overvoltage category:	II
Degree of pollution:	2
Ingress protection code:	IP43 (balances with readability of 0.01 g or higher, excluding portable balances)

#### Note

Stated IP is only achieved when the balance is ready for operation. The protective cover must be installed, and the caps must cover the interface connections.

Standards for safety and EMC:	See Declaration of Conformity
Range of application:	Use only indoors in dry locations

## Environmental conditions

The limit values apply when the balance is used under the following environmental conditions:

Height above mean sea level:	Up to 5000 m
Ambient temperature:	+10 – +30 °C (excluding large balances) +5 – +40 °C (large balances)
Temperature change, max.:	5 °C/h
Relative humidity:	30 – 70%, non-condensing
Acclimatization time:	Recommendation: Up to <b>4 hours</b> for precision balances, or up to <b>8 hours</b> for analytical balances. These values apply after placing the balance in the same location where it will be put into operation.

### Note

The acclimatization time depends on the readability of the balance, and on the environmental conditions.

Warm-up time:	At least <b>30 minutes</b> for precision balances, or <b>60 minutes</b> for analytical balances. These values apply after connecting the balance to the power supply. When switched on from standby, the balance is ready for operation immediately.
---------------	--

The balance can be used under the following environmental conditions. However, the weighing performances of the balance may be outside the limit values:

Ambient temperature:	+5 °C – +40 °C
Relative humidity:	20% to max. 80% at 31 °C, decreasing linearly to 50% at 40 °C, non-condensing

The balance can be disconnected and stored in its packaging under the following conditions:

Ambient temperature:	-25 – +70 °C
Relative humidity:	10 – 90%, non-condensing

## 7 Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties, the content of this regulation must also be related.

## 8 Compliance Information

National approval documents, e.g., the FCC Supplier Declaration of Conformity, are available online and/or included in the packaging.

► [www.mt.com/ComplianceSearch](http://www.mt.com/ComplianceSearch)

For further information, consult the Reference Manual (RM).

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► [www.mt.com/contact](http://www.mt.com/contact)

## 1.3 Akronymmer og forkortelser

Originalt udtryk	Oversat udtryk	Forklaring
ABS/PC		Acrylonitrile butadiene styrene / polycarbonate (polymer blend)
AC		Alternating Current
ASTM		American Society for Testing and Materials
DC		Direct Current
EMC		Electromagnetic Compatibility
FCC		Federal Communications Commission
GWP		Good Weighing Practice
ID		Identification
IP		Ingress Protection
LPS		Limited Power Source
MT-SICS		METTLER TOLEDO Standard Interface Command Set
NA		Not Applicable

OIML	Organisation Internationale de Métrologie Légale (International Organization of Legal Metrology)
PA 12	Polyamid 12
PBT	Polybutylene terephthalate
PET	Polyethylene terephthalate
POM	Polyoxymethylene
RM	Reference Manual
SOP	Standard Operating Procedure
TDNR	Type Definition Number
TPE	Thermoplastic elastomer
UM	User Manual (Brugervejledning)
USB	Universal Serial Bus
USP	United States Pharmacopeia

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© 07/2024 METTLER TOLEDO. All rights reserved.