

# NewClassic Balances ME Models

## Simple Weighing

**WEIGH** → WEIGH Activate simple weighing  
• Exit application  
0.00 g

**Tare**  
0.00 g  
→0/T← → 12.01 g  
0.00 g NET

**Switching Units**  
22.00 g  
\* 22.00 g<sup>M</sup> Recall  
• if activated  
0.78 oz Unit 2  
• if activated  
• if different from Unit 1

## Weighing Applications

**F** → COUNT

**APPL LIST** → PERCENT %

CHECK

STAT

**←** ← FORMULA

TOTAL

DYNAMIC

FACTOR.M Fx

FACTOR.J F+

DENSITY

## Adjustments

**Cal** → External Adjustment Procedure

ADJUST →

← ← ADJUST

← ← 1000.00 g

← ← 1000.00 g

← ← 0.00 g

← ← ADJ.DONE

← ← 0.00 g

## Operation Keys

Indicates active application →

Weighing-in aid →

Indicates unstable values →

1 C ΔΔ

2 F

3 →0/T←

4 Cal

5 Menu

Key Functions			
No.	Key	Press briefly (less than 1.5 s)	Press and hold (longer than 1.5 s)
1	C ΔΔ	<ul style="list-style-type: none"> <li>Cancel or leave menu without saving</li> <li>One step back in the menu</li> </ul>	<ul style="list-style-type: none"> <li>Select the simple weighing application</li> <li>Exit application</li> </ul>
2	F	<ul style="list-style-type: none"> <li>Printout display value</li> <li>Transmit data</li> <li>To navigate backwards in the menu and menu selection</li> <li>Decrease parameters in menu and applications</li> </ul>	<ul style="list-style-type: none"> <li>Open the application list for selecting an application</li> </ul>
3	→0/T← ⏻	<ul style="list-style-type: none"> <li>Zero/Tare</li> <li>Switch on</li> </ul>	<ul style="list-style-type: none"> <li>Switch off into standby mode</li> </ul>
4	↻ Cal	<ul style="list-style-type: none"> <li>With entries, scroll down</li> <li>To navigate forward menu topics or menu selections</li> <li>To toggle between unit 1, recall value (if selected), unit 2 (if different from unit 1) and the application unit (if any)</li> <li>Increase parameters in menu or applications</li> </ul>	<ul style="list-style-type: none"> <li>Select adjustment (calibration)                             <ul style="list-style-type: none"> <li>– with internal weight *</li> <li>– with external weight</li> <li>– customer fine adjustment *</li> </ul> </li> <li>* On models with internal weight only</li> </ul>
5	↵ Menu	<ul style="list-style-type: none"> <li>Enter or leave menu selection</li> <li>To enter application parameter digit and switch to next parameter digit</li> <li>To accept parameter in menu selection</li> </ul>	<ul style="list-style-type: none"> <li>Enter or leave menu (parameter settings)</li> <li>To store parameter</li> <li>To accept numeric input in applications</li> </ul>

# Menu Map

## Main Menu

BASIC	←	Basic Menu
ADVANCE.	←	Advanced Menu
INT.FACE	←	Interface Menu
PROTECT	←	Protection

## Protection

Menu	Selection
PROTECT	← OFF
	← ON

## Basic Menu

Topic	Selection
DATE	← 01.01.11
TIME	← +1H
	← -1H
	← SET.TIME 12:00
UNIT 1	← OFF
	← ON
UNIT 1	← g
	← kg
UNIT 2	← g
	← mg
SET ID	←
PRT.MENU	←
RESET	← NO?
	← YES?

## Advanced Menu

Topic	Selection
ENVIRON.	← STB.
	← UNSTAB.
	← STABLE
ADJ.LOCK	← OFF
	← ON
DATE.FRM	← 00MMYY
	← MM/00YY
	← Y-MM-00
	← 0MMYY
	← MMYY
TIME.FRM	← 24MM
	← 12MM
	← 24MM
	← 12MM
RECALL	← OFF
	← ON
STANDBY	← A.ON
	← A.OFF
	← 60
LIGHT	← 3L.ON
	← 3L.OFF
A.ZERO	← ON
	← OFF
ZERO.RNG	← 21.00 g
SRV.ICON	← ON
	← OFF
SRV.RST	← NO?
	← YES?

## Interface Menu

Topic	Selection
RS232	← PRINTER
	← PRT.STAB
	← PRT.AUTO
	← PRT.ALL
PC.DIR	← PRT.STAB
	← PRT.AUTO
	← PRT.ALL
HOST	← SND.OFF
	← SND.STB
	← SND.CONT
	← SND.AUTO
	← SND.ALL

Topic	Selection
HEADER	← 2DISP
	← NO
	← PARTIM
	← PARTBL
SINGLE	← NET
	← S/T/N
SIGNL	← OFF
	← ON
LN.FEED	← 00
ZERO.PRT	← OFF
	← ON
COMSET	← MT-SICS
	← SART
AUD	← 9600
	← 4800
	← 600...38400
BIT.PAR.	← 8/NO
	← 7/NO
	← 7/MARK
	← 7/SPACE
	← 7/EVEN
	← 7/ODD
STOPBIT	← 1 BIT
	← 2 BITS
HD.SHK	← %ON.%OFF
	← RTS.CTS
	← OFF
RS.TX.EOL	← CR LF
	← CR
	← LF
	← TAB
RS.CHAR	← 13M.DOS
	← ANSL.WIN
INTERVL	← 00000

### Legend

- Factory setting
- Only those weight units allowed by the appropriate national legislation are selectable
- Not available with approved models
- Not available with approved models with e = d
- Settings are automatically set for the 2<sup>nd</sup> display
- Only visible if "PRINTER" is selected
- Only visible if "HOST" is selected
- Only visible if "PC-DIR" is selected
- Only visible if "PRT.AUTO" is selected

[www.mt.com/newclassic](http://www.mt.com/newclassic)

For more information

Mettler-Toledo AG  
Laboratory & Weighing Technologies  
CH-8606 Greifensee, Switzerland

Subject to technical changes  
© Mettler-Toledo AG 07/2012  
30045679A en



\* 3 0 0 4 5 6 7 9 \*