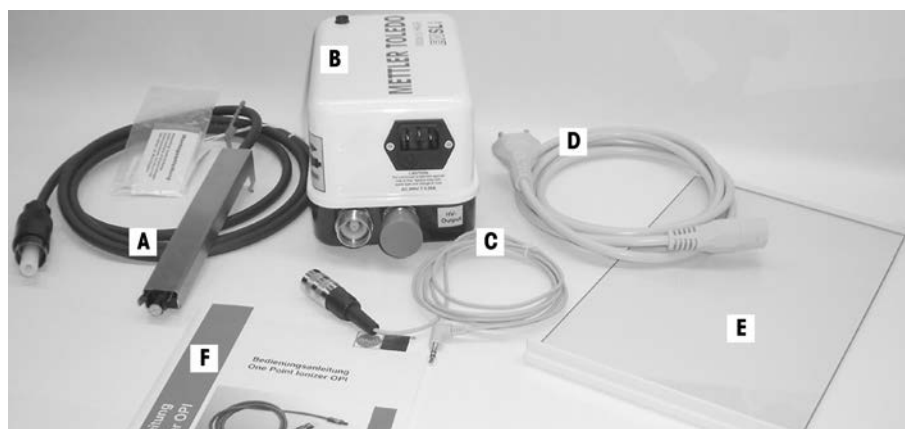
Installation Instructions **Antistatic Kit for XP Analytical Balances**  
 Montageanleitung **Antistatik Kit für XP Analysenwaagen**  
 Notice de montage **Kit antistatique pour les balances d'analyse XP**  
 Instrucciones de montaje **Kit antiestático para balanzas analíticas XP**  
 Istruzioni di montaggio **Kit antistatico per bilance analitiche XP**

**Components**

- |                               |                                 |
|-------------------------------|---------------------------------|
| <b>A</b> Single-point ionizer | <b>B</b> AC adapter             |
| <b>C</b> Control cable        | <b>D</b> Power supply cable     |
| <b>E</b> Side door            | <b>F</b> Operating Instructions |

**Note**

Position the single-point ionizer on the left or right side of the balance.  
The AC adapter can have two single-point ionizers connected to it.

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### CAUTION

#### Damage of Device

The power cable used to connect the ionizer must have a ground wire and be plugged into an outlet with a functioning ground connection.

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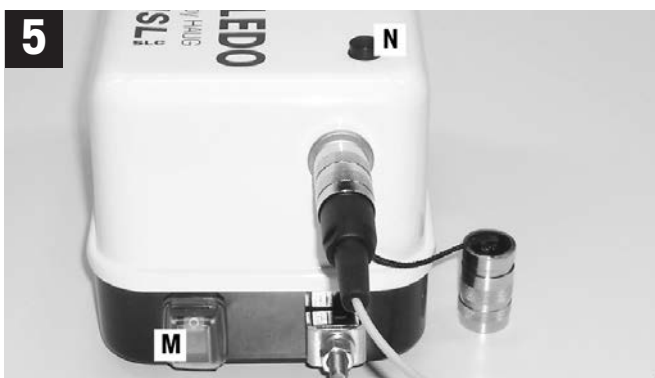
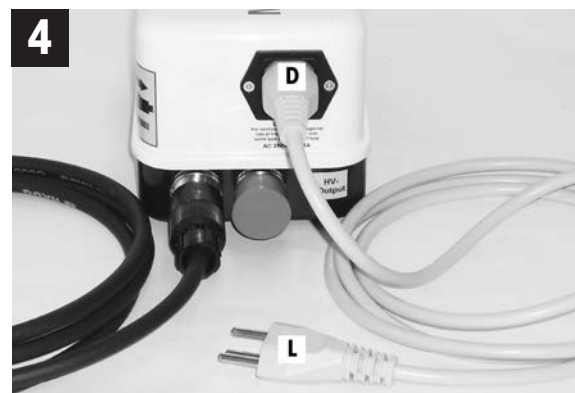
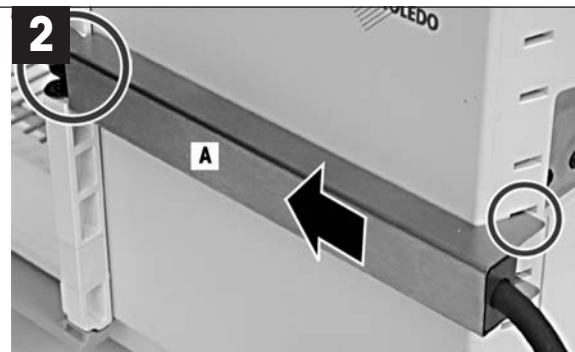
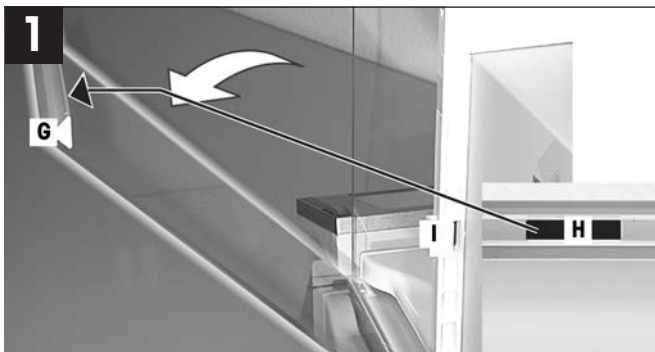
## English

- Swap the side door (G) for the door (E) marked with "Conducting Glass" (H) supplied with the antistatic kit.
  - Pull the middle clip (I) (recommended position) off.
- Push the point ionizer (A) in from the back. If necessary, slightly bend the strips so that the ionizer is held firmly.
- Screw the ionizer cable (J) tightly onto the AC adapter.
  - Screw in the control cable (C).
  - Plug the control cable into Aux 1 or Aux 2 (K).
- Plug in power cable (D).
  - Connect power cable (L) to power supply.
- Switch On/Off (M).
  - ⇒ Control lamp (N) lit when the ionizer is active.

Information about settings for the antistatic kit is contained in the operating instructions for XP analytical balances.

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[www.mt.com/excellence](http://www.mt.com/excellence)

For more information

**Mettler-Toledo AG**

Laboratory & Weighing Technologies  
 CH-8606 Greifensee, Switzerland  
 Phone +41 44 944 22 11  
 Fax +41 44 944 30 60

Subject to technical changes  
 © Mettler-Toledo AG 09/2020  
 30038450B en, de, fr, es, it