M80 SM **Sensor Mount Transmitter**

Instruction Manual

cs	Návod k obsluze	2
da	Brugsanvisning	8
de	Betriebsanleitung	14
en	Instruction manual	22
es	Manual de instrucciones	27
fi	Käyttöohje	34
fr	Instructions d'utilisation	40
hu	Használati utasítás	47
it	Istruzioni per l'uso	54
ja	取扱説明書	61
ko	사용 매뉴얼	67
nl	Gebruikershandleiding	73
pl	Instrukcja obsługi	80
pt	Manual de instruções	87
ru	Инструкция по эксплуатации	94
sv	Brugsanvisning	102
th	คู่มือคำแนะนำ	108
zh	说明书	115

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1 Regulatory Approval

The measuring system is in conformity with the statutory requirements of the EC Directives. METTLER TOLEDO confirms successful testing of the device by affixing to it the CE mark.

2 Environmental Protection

Waste electrical products should not be disposed of with household waste. Please recycle where facilities exist. Check with your Local Authority or retailer for recycling advice.

3 Safety Instructions

3.1. Definition of Equipment and Documentation Symbols and Designations

WARNING: POTENTIAL FOR PERSONAL INJURY

CAUTION: Possible instrument damage or malfunction

NOTE: Important operating information

On the transmitter or in the manual text indicates: Caution and/or other possible hazard including risk of electric shock (refer to accompanying documents).

3.2. General safety instructions and warnings

 **NOTE: Be sure to read and observe the following instructions.**

The following is a list of general safety instructions and warnings. Failure to adhere to these instructions can result in damage to the equipment and/or personal injury to the operator.

- The M80SM transmitter should be installed and operated only by personnel familiar with the transmitter and who are qualified for such work.
- The M80SM transmitter must only be operated under the specified operating conditions.
- Repair of the M80SM transmitter must be performed by authorized, trained personnel only.
- With the exception of cleaning procedures, as described in this manual, the M80SM transmitter must not be tampered with or altered in any manner.
- METTLER TOLEDO accepts no responsibility for damage caused by unauthorized modifications to the transmitter.
- Follow all warnings, cautions, and instructions indicated on and supplied with this product.
- Install equipment as specified in this instruction manual. Follow appropriate local and national codes.
- If this equipment is used in a manner not specified by the manufacturer, the protection provided by it against hazards may be impaired.

 **NOTE: Important operating information**
Before connecting the device to a supply unit, make sure that its output voltage cannot exceed 30 VDC. Do not use alternating current or main power supply!

3.3. Warnings

- Installation of cable connections and servicing of this product require access to shock hazard voltage levels.
- Electrical installation must be in accordance with the National Electrical Code and/or any other applicable national or local codes.

 **NOTE: Process Upsets:** Because process and safety conditions may depend on consistent operation of this transmitter, provide appropriate means to maintain operation during sensor cleaning, replacement or sensor or instrument calibration.

 **WARNING: Do not install the M80SM transmitter in hazardous locations!**

3.4. Correct disposal of the unit

When the transmitter is finally removed from service, observe all local environmental regulations for proper disposal.

4 Introduction

Thank you for buying this METTLER TOLEDO equipment. The M80SM is a direct mount transmitter for METTLER TOLEDO ISM[®] (Intelligent Sensor Management) 1-wire sensors for digital integration into benchtop biocontrollers via an RS485 MODBUS RTU interface. For the support of MODBUS RTU programming please download the M80SM MODBUS Programmers Guide from the documentation section on www.mt.com/M80.

If configuration of an M80SM transmitter is necessary use the METTLER TOLEDO M80 Transmitter Configuration Tool (TCT) software and the METTLER TOLEDO Transmitter Configuration Box (see chapter 9 for ordering information). Please download the TCT software from the documentation section on www.mt.com/M80.

The M80SM stores installation point-specific MODBUS parameters. So there's no need to re-configure the transmitter when a 1-wire sensor is replaced.

5 Technical data M80 SM

Measurement parameters	1-wire sensors: pH, amperometric Dissolved Oxygen, Dissolved CO ₂ , Conductivity
ISM	Full ISM functionality
Power supply	24 VDC (min. 100 mA) / 8 – 30 VDC (min. 2 W)
Operating temperature	–15 to +60 °C
Relative humidity	5 ... 95 % rH (non-condensing)
Protection class	IP65

6 Connections

The M80SM 1-wire has one connector to fit any METTLER TOLEDO 1-wire ISM sensor (AK9 connector).

NOTE: Important operating information
The M80SM transmitter cannot be auto-claved! Disconnect the M80SM transmitter from the sensor before autoclaving the sensor!

The 5-pin M12 connector is connected to a combined 5-pin data/24 VDC power supply cable that is available from METTLER TOLEDO in different lengths – see chapter 9 for ordering information.

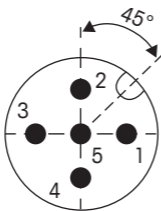
6.1. Connection to controller

 **NOTE:** For connecting the M80SM to a controller MODBUS RTU interface and a 24 VDC power supply use a METTLER TOLEDO standard M12 5-pin data cable with open ends.

Pin and cable assignment M12 connector (male) and 5-pin data cable:

Male M12

connector pin	Open cable end	Function
1	brown	24 VDC+
4	black	24 VDC-
3	blue	RS485- (B)
2	white	RS485+ (A)
5	gray	ground
	yellow	shield

Male M12 connector on M80SM transmitter

7 MODBUS configuration

MODBUS specific settings (baud rate, device address, etc.) can be adjusted when the M80SM is connected to the Transmitter Configuration Tool PC Software via the Transmitter Configuration Box with USB adapter. Please refer to the Transmitter Configuration Box manual for details.

Please download the M80 TCT software from www.mt.com/M80.

8 Cleaning/storage

Clean the surfaces of the M80SM transmitter with a soft damp cloth, then dry the surface with a cloth, carefully. Do not use any solvents or water to clean connecting parts of connectors!

For storage, the device should be clean and dry.

9 Ordering information

Item	Ordering number
M80SM Transmitter	30 530 566
Transmitter Configuration Box	30 530 567
5-pin data cable 2 m (6.6 ft)	52 300 379
5-pin data cable 5 m (16.4 ft)	52 300 380
5-pin data cable 10 m (32.8 ft)	52 300 381

ISM is a registered trademark of the METTLER TOLEDO Group in Switzerland, Brazil, USA, China, European Union, South Korea, Russia and Singapore.

All other trademarks are the property of their respective holders.

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All rights reserved.